	TECHNICAL DELIVERY CONDITION 50145670	Page: 1 of 2 Date: 02.12.2020
	Article name: mySPIN6 Instruction Manual	Internal note: mySPIN6 Instruction Manual.pdf
		Editor:
		Approval:

Article name: MYSPIN 6 INSTRUCTION MANUAL

Editor: Gayathri Duggirala

Date: 02.12.2020

Place: R&D IEC HYD

Table of Contents

1	Description	2
2	Specifications	2
3	Attachment	2

History

f	TBD	Correction of product name, re-translation from EN	02.12.2020	Gayathri Duggirala
e	219460	Updated as per ECR 219460	26.08.2020	Gayathri Duggirala

Rev.	ECR No.	Description	Date	Name CHKD
f	TBD	Correction of product name, re-translation from EN	12.02.2020	Gayathri Duggirala

	TECHNICAL DELIVERY CONDITION 50145670	Page: 2 of 2 Date: 02.12.2020	
	Article name: mySPIN6 Instruction Manual	Internal note: mySPIN6 Instruction Manual.pdf	
			Editor:
			Approval:
Gayathri Duggirala			

1	Description
	Refers to the instruction manual for the mySPIN 6 in English, French, Italian, German and Spanish.
2	Specifications
	Only relevant, if printed! <ul style="list-style-type: none"> • Print in upright format. Custom page size: 4.9 inch x 8.50 inch. • 2 pages cover • Cover bright white 160 g/m² • Content bright white 80 g/m² • Double-sided print, colored • Stapled
3	Attachment
	User Instruction sheet in PDF format.

Rev.	ECR No.	Description	Date	Name CHKD
f	TBD	Correction of product name, re-translation from EN	12.02.2020	Gayathri Duggirala

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ThermoFisher
SCIENTIFIC

Contents

Safety Precautions	1
Specifications	5
Environmental Ratings.....	5
Technical Features	5
Electrical	5
Accessories	6
Operation.....	8
Operating the mySPIN 6 in a Lab	8
Operating the mySPIN 6 in a Remote	10
Environment.....	10
Rotor Replacement	10
Maintenance	11
Troubleshooting	12
Disposal	12
Regulatory Information	13
Warranty	14
Contact Information	15

Safety Precautions

The Thermo Scientific™ mySPIN 6 is a personal mini centrifuge designed for laboratories in the bioscience, medical, and chemistry fields. It may be used with biohazardous and/or hazardous materials. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instrument's suitability to the task.

Before starting the centrifuge for the first time, read this instruction manual carefully. This manual contains important information, safeguards, and operating instructions.

CAUTION Never use this product in any manner inconsistent with these instructions.

CAUTION This product is intended for separating aqueous solutions in approved test tubes compatible to the RCF of this unit.

CAUTION This product is not intended to centrifuge very dense materials. Do not use materials with a density greater than 1.2 g/cm³ (1.2 g/mL).

CAUTION Use only the power adapter delivered with the product. Use of other power adapters will void warranty.

CAUTION Do not put hands in the rotor area unless the rotor is completely stopped.

CAUTION Only use the originally supplied rotors, AC adapters, tube adapters and any other spare parts delivered with the unit. These parts have been designed to safely meet the maximum RCF that the centrifuge is able to deliver. **DO NOT USE ROTORS OR COMPONENTS FROM ANY OTHER MANUFACTURER.**

CAUTION Always make sure the rotors are securely seated.

CAUTION Observe general laboratory safety precautions and regulations.

CAUTION Do not attempt to operate with the cover removed or open.

CAUTION The rotor and lid must always be securely fastened when in operation.

CAUTION Do not use damaged rotors.

CAUTION Do not move unit while the rotor is spinning.

CAUTION Do not attempt to stop the rotor while the unit is running. Doing so may cause injury or the unit to fail and will void the warranty.

CAUTION mySPIN 6 mini centrifuge shall not be operated in wet conditions, including but not limited to rain, snow, immersion in water. Injury due to shock or fire may result.

CAUTION The rotor must be loaded symmetrically. Operating this unit with an unbalanced rotor will cause severe vibration and damage, which is not covered by warranty.

CAUTION Do not fill tubes while inserted in rotor. Spilled liquid may harm unit.

CAUTION Close tube lids before starting centrifuge. Open lids can be ripped off during centrifugation and damage the unit. Open lids can cause fluid to be dispersed inside the centrifuge.

CAUTION Do not insert non-standard tubes or other unfamiliar articles into the rotor.

CAUTION Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this instrument.

CAUTION Do not use solvent or flammable liquids near this or other electrical equipment.

CAUTION Do not centrifuge materials that in combination or alone are flammable, explosive, or corrosive, or may cause any other hazard.

CAUTION Do not operate in a hazardous or flammable environment.

CAUTION Do not use any solvent on the unit that may attack plastic or rubber or cause cracks in the rotor.

CAUTION Ensure the rotor is protected from corrosion and mechanical damage. The rotor must be cleaned with a pH-neutral cleaning liquid.

CAUTION Before using any cleaning or decontamination methods except those recommended in this manual, contact Thermo Fisher Scientific to ensure that the proposed method will not damage the device.

CAUTION If a single level of hazard protection is required, make sure to use containers that are tested to Risk Group II or above and that more than one level of protection is provided in the case of materials of a higher Risk Group.

CAUTION Should the unit fail to work to specification, stop centrifuging immediately and allow the rotor to come to a complete stop. Clean and troubleshoot the product according to the instructions under “**Troubleshooting**” before any further use of the product. Contact an authorized dealer or the manufacturer if trouble persists.

CAUTION Always work in a manner which endangers neither the user nor any other person.

CAUTION Repairs are to be performed by trained and authorized service personnel only. Contact Thermo Fisher Scientific.

CAUTION Opening the instrument housing or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Thermo Fisher Scientific for warranty service.

CAUTION Do not use units that have not been correctly installed or repaired.

CAUTION Do not attempt to disassemble or modify this product.

CAUTION Store at room temperature in a dry area. Do not expose to sunlight, moisture, or extreme temperatures for prolonged periods of time.

WARNING The laboratory procedure shall ensure that no person or hazardous substances are present within a 12" (30 cm) zone around the centrifuge when operating.

WARNING If this product is not used as recommended by the manufacturer, the overall safety will be impaired.

WARNING Check MSDS, wear required Personal Protective Equipment, and observe all applicable local and national regulations before dispensing and disposing of hazardous materials.

WARNING This product does not contain bio-seals as per IEC/EN/CSA 61010-2-020 and cannot provide any level of containment in case of a spill or release of toxic, radioactive, or pathogenic micro-organisms. Thus these materials are not recommended to be used with this product.

Specifications

Environmental Ratings

Altitude:	up to 2000 m
Operating Temperature:	0 °C to 40 °C (32 °F to 104 °F)
Storage Temperature:	-20 °C to 40 °C (-4 °F to 104 °F)
Maximum Relative Humidity:	80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40 °C
Relative Humidity during operation:	10% to 90% (non-condensing)
Relative humidity during storage:	10% to 90%

Technical Features

Capacity:	Standard rotor 6 x 2.0 mL PCR strip rotor 2 x 8 x 0.2 mL
Maximum speed:	6,000 rpm
Maximum RCF:	2,000 x g
Dimensions:	153 mm x 128 mm x 104 mm (L x W x H)
Weight:	1.6 lb (0.73 kg) (with rotor) 1.9 lb (0.86 kg) (with power adapter)

Electrical

mySPIN 6

Input: 12 V DC

Current (under load): Approx. 0.6 A

Power: Approx. 7.3 W

Power Adapter rating

Input: 100–240 V AC 50/60 Hz, 1.0 A

Output: 12 V DC, 2.2 A

12 V Car Adapter

Input / Output: 12–13.8 V

Rating: 12 V DC

Accessories

Cat. No	Description
75004061	mySPIN 6 mini centrifuge Includes: <ul style="list-style-type: none">• 6 x 1.5 / 2.0 mL tube rotor• 16 x 0.2 mL tube rotor (singles or 2 x 8 strip)• 0.2 mL tube adapters (set of 6)• 0.5 mL tube adapters (set of 6)• Storage case for rotors and adapters• Tube rack• Universal power adapter 100–240 V 50 / 60 Hz, with 4 inter-changeable plug types
75004085	mySPIN 6 Portable Kit Complete
Replacement Parts	
75004062	6 x 1.5 / 2 mL circular rotor
75004063	16 x 0.2 mL tube rotor (singles or 2 x 8 strip)
20290729	Replacement O-ring
50146145	Tube adapter inserts for 0.2 / 0.5 mL (12 per pack; 6 of each size)
75004064	Power adapter and universal power cord set
75004086	mySPIN 6 Transport Box with inserts
75004087	mySPIN 6 12 V Car Adapter

mySPIN 6 Portable Kit (75004085):

The mySPIN 6 Portable Kit is intended for separating samples outside the laboratories or in a remote location. The kit includes:

- mySPIN 6 mini centrifuge complete with instruction manual
- Clear lockable transport box
- Set of dividers with integral tube rack, dividers designed to fit mySPIN 6
- 12 V automobile power adapter, for connection from cigarette lighter socket to mySPIN 6 mini centrifuge.

mySPIN 6 Transport Box with Inserts (75004086):

The transport box is intended to accommodate the mySPIN 6 mini centrifuge with other necessary accessories, like rotor and tube rack to carry them out from the laboratories for field work.

- Dimensions: 380 mm x 196 mm x 160 mm (L x W x H)
- Weight: 1.2 kg
- Autoclavable
- Watertight (leak-proof), silicone seal
- Leak proof when submerged in 1 meter of water
- Supports up to 4 kg of weight

The 6 inserts inside the transport box are shown in the figure below.

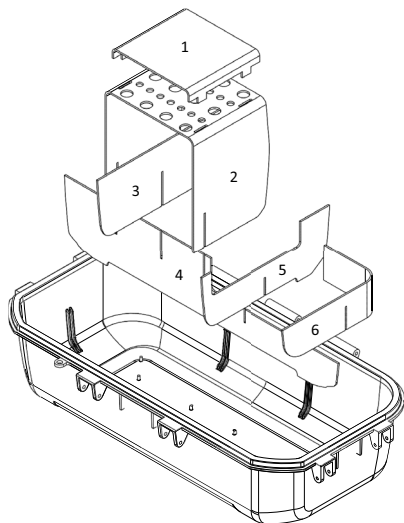
Figure 1. Inserts of Transport Box

You may add or remove the inserts as per your requirements.

12 V Car Adapter (75004087):

The 12 V car adapter is used to power up the mySPIN 6 unit and perform the sample separation in the field.

- Cable length: 1,500 mm
- 24 awg / 2 C cable
- Plug fit in different cars with a cigarette lighter socket

Operation

Operating the mySPIN 6 in a Lab

1. Remove the contents from the package, examining them carefully for breakage, defects, or missing parts. Save the packaging.

Notice The unit shall be discarded if the locks or housing are broken or cracked. Follow your facility or region policies for the disposal of biological hazardous materials.

2. Place the mySPIN 6 on a sturdy, level surface away from the edge of the counter and other moving equipment. Make sure the power cord is clear of hot surfaces and other hazards.

WARNING Spin balanced loads only. Do not place the tubes asymmetrically nor load unequal volume tubes. Improper placement of tubes will lead to insufficient centrifugation and may cause injury or accident.

3. Prior to every use, inspect unit for cracks, missing pieces, and/or any damage that would result in an impaired performance and notify Thermo Fisher Scientific BEFORE use.
4. Verify that the power switch is in the OFF position. Plug the power adapter cord into the rear of the unit and the power cord into an approved outlet. Make sure that the power cord connector can be readily grasped at the rear of the device so that it can be removed to easily disconnect it from the centrifuge.
5. Turn power switch ON. If there is a smooth whirring sound and the unit accelerates with little or no vibration, mySPIN 6 is ready to use. If there are loud or unusual sounds, or excessive vibration, DO NOT OPERATE. Contact Thermo Fisher Scientific Technical Support.
6. On benchtop prepare samples tubes in rack so that tubes are filled to equal levels. Close tube lids.
7. To open lid, firmly tilt lid toward rear of unit.

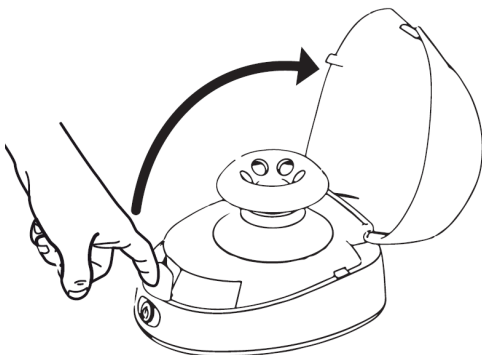
Figure 2. Open lid of mini centrifuge

WARNING The laboratory procedure shall ensure that no person or hazardous substances are present within a 12" (30 cm) zone around the centrifuge when operating.

- Tubes of equal weight and size shall be placed opposite each other. Use sample or water in tube to provide balance. Proper sample balancing illustrated below.
- Do not force tubes into the rotor. Use adapters and correctly sized rotors for your application.

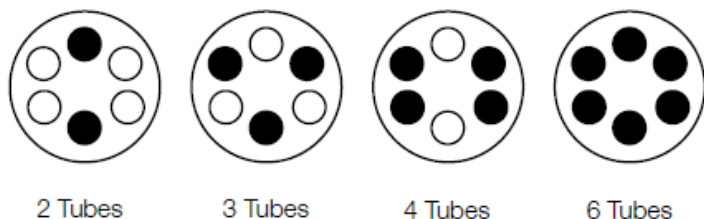
Figure 3. SnapSpin Circular Rotor

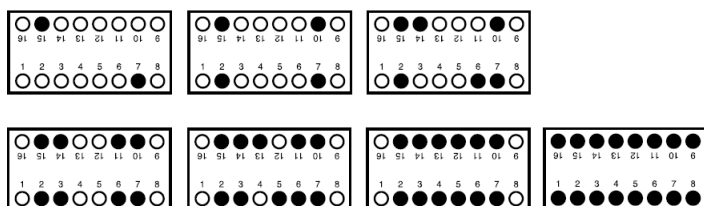
Figure 4. SnapSpin Strip Tube

- Close lid securely. A properly closed lid will grasp the front silicon tab keeping the lid closed. If a clicking sound is heard while the unit is operating, the lid is not fully closed. Do not attempt to operate with the lid removed or open. Turn power switch ON. Unit will not operate with lid open. If there is a smooth whirring sound and the unit accelerates with little or no vibration, mySPIN 6 is ready to use. If there are loud or unusual sounds, or excessive vibration, DO NOT OPERATE. Contact the Thermo Fisher Scientific Technical Support.
- To stop centrifugation, press power switch and wait until the rotor has completely stopped rotating before opening the lid.

WARNING Do not open lid until rotor has completely stopped! Avoid severe personal injury or property damage from moving parts.

Operating the mySPIN 6 in a Remote Environment

1. Place the mySPIN 6 on a sturdy, level surface away from the edge of the counter and other moving equipment.
2. Verify that the power switch is in the OFF position.
3. Plug the 12 V car adapter into the car cigarette lighter socket.
4. Press the power button  on the centrifuge.
5. Follow step 7 to step 10 from the procedure above in **Operating the mySPIN 6 in a Lab**.
6. Remove the adapter from the car.

CAUTION Prevent any twists or bends to the adapter cable to avoid any damages.


Notice The unit must be sealed securely and locked when carried with samples enclosed.

Rotor Replacement

To remove SnapSpin™ rotor: Press center button firmly and at the same time lift rotor upwards.

To install SnapSpin™ rotor: Set rotor onto the hex shaft. Press center button firmly and downward. Rotor shall latch with a click sound.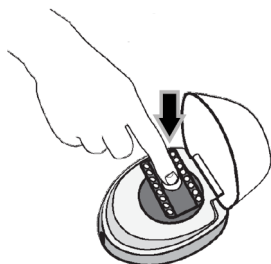

Maintenance

mySPIN 6 is normally maintenance-free. Clean unit only when it is not plugged into an electrical outlet. When necessary, the housing and rotor can be wiped using a damp cloth and a mild, non-corrosive detergent.

WARNING Do not use any solvent on the unit that may attack plastic or cause cracks in the rotor or the unit. Avoid exposure to UV radiation, either natural or man-made. In addition to cracks; discoloration, softening, or crazing of the plastic materials can occur. Corrosion can also occur to the metal parts which are part of the SnapSpin™ rotor feature. This damage can weaken them. Ensure the rotor is protected from corrosion and mechanical damage. The rotor must be cleaned with a neutral cleaning liquid.

Notice Ensure all parts are dry prior to operation.

The transport box can be cleaned using soap, water, bleach mixture, IPA & Ethanol. It is autoclavable at 121°C for 15 minutes.

WARNING Higher temperatures or autoclaving duration may damage the transport box. The box shall be unlocked and open during autoclaving with the silicone seal left in place.

Troubleshooting

Table 1. Troubleshooting

Problem	Possible Cause	Solution
Not working No light	Loose connection of power cord / adapter	Check power cord / adapter
	Front ON / OFF switch	Press front ON / OFF switch
	Lid not closed	Close Lid
Excessive vibration	Rotor not installed correctly	Reinstall rotor
	Rotor not balanced correctly	Check for tubes of equal weight and location (see Operation)
	Soft or uneven surface	Place the mySPIN 6 on a sturdy and level surface.
Noise clicking sound	Lid not fully closed	Close Lid engaging silicon tab with cover
Rotor falls off	Rotor not attached	Reattach rotor (See Rotor Replacement)

Disposal

The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.

If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact us for more information on the environmental performance of our products.

Regulatory Information

This mark is the confirmation that the unit conforms to the EU requirements.

Conforms to:

ANSI/UL Std. 61010-1

IEC/EN 61010-1

IEC/EN 61010-2-020

CAN/CSA C22.2 No. 61010-1.

Warranty

IN NO EVENT WILL THERMO FISHER SCIENTIFIC OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

Limited Warranty

Thermo Fisher Scientific warrants that mySPIN 6 will be free from defects in workmanship and material for 3 years from the date of purchase.

Thermo Fisher Scientific provides one year warranty for the transport box and 12 V car adapter from date of purchase.

If you believe that there is a defect in the product, you must, during the warranty period, notify Thermo Fisher Scientific, provide proof of purchase, and return the product to Thermo Fisher Scientific with a Return Authorization form.

If Thermo Fisher Scientific is properly notified and, after inspection, confirms that there is a defect and the warranty period has not expired, Thermo Fisher Scientific will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, THERMO FISHER SCIENTIFIC MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. THERMO FISHER SCIENTIFIC WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

ACCESSORY PARTS, SUCH AS RUBBER PARTS, THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (mySPIN 6 IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)

ANY PRODUCT THAT HAS BEEN, IN THERMO FISHER SCIENTIFIC SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN THERMO FISHER SCIENTIFIC.

ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.

Register your warranty online at

www.thermoscientific.com/warranty.

thermo scientific

Contact Information

Thermo Fisher Scientific products are backed by a global technical support team ready to support your applications. Visit www.thermofisher.com or call:

Countries	Sales	Services
North America	+1 866 984 3766	(800) 438-4851
India	1800 22 8374, +91 22 6716 2200	+91 22 6716 2200
China	+800 810 5118, +400 650 5118	+8621 68654588
Japan	+81 3 5826 1616	+81 3 3816 3355
Australia	+61 39757 4300	1 300 735 292
Austria	+43 1 801 40 0	+43 1 801 40 0
Belgium	+32 53 73 42 41	+32 2 482 30 30
France	+33 2 2803 2180	+33 2 2803 2180
Germany	0800 1 536 376, +49 6184 90 6000	0800 1 536 376
Italy	+32 02 95059 552	+39 02 95059 552, 432 254 375
Netherlands	+31 76 579 55 55	+31 76 571 4440
Nordic/Baltic/CIS	+358 9 329 10200	+358 9 329 100
Russia	+7 812 703 4215	+7 812 703 4215
Spain/Portugal	+34 93 223 09 18	+34 93 223 09 18
Switzerland	+41 44 454 12 22	+41 44 454 12 12
UK/Ireland	+44 870 609 9203	+44 870 609 9203
New Zealand	+64 9 980 6700	+64 9 980 6700
Other Asian Countries	+852 2885 4613	+852 2885 4613
Countries not listed	+49 6184 90 6000	+49 6184 90 6000

Distributor: Thermo Electron LED GmbH
Zweigniederlassung Osterode
Am Kalkberg, 37520 Osterode am Harz
Germany

Find out more at thermofisher.com/Centrifuge

© 2016–2020 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. Heathrow Scientific is a registered trademark of Heathrow Scientific. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Consult your local sales representative for details. Shown pictures within the manual are examples and may differ considering the set parameters and language.

50145670 (English) is the original instruction manual.

Manufacturer: Heathrow Scientific
620 Lakeview Parkway, Vernon Hills
IL, 60061
USA

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EN NINGÚN CASO LA OBLIGACIÓN DE THERMO FISHER SCIENTIFIC LLC SEGÚN ESTA GARANTÍA SUPERARÁ EL PRECIO DEL PRODUCTO.
Distribuidor: Thermo Electron LED GmbH