MB90 / MB120 Moisture Analyzer User Guide

EN

ES

FR

DE

IT

PT

SV

DA

PL

NL

CZ

HU

1. SAFETY INFORMATION

Definition of Signal Warnings and Symbols

WARNING	For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
CAUTION	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
Attention	For important information about the product.
Note	For useful information about the product.

Warning Symbols

General hazard

Electrical shock

Hot surface

Inflammable or explosive substance

Toxic substance

Acid / Corrosion

Safety Precautions

CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

Intended Use

Use the instrument exclusively for <weighing/moisture determination/etc.> as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

The Moisture Analyzer works with heat!

- Never place flammable materials on, below or next to the Analyzer.
- Use caution when removing a test sample. The sample, the sample chamber, the heating element and the surrounding areas may be very hot and can cause burns.

Some samples require special care!

- Should there be any uncertainty regarding the safety of a substance, perform a careful risk analysis. In such cases, never leave the Analyzer unattended.

- Fire or explosion: Substances which contain solvents or release flammable or explosive vapors when heated. With such samples, work at drying temperatures low enough to prevent the formation of flames or an explosion.

- Poisoning or burning: Substances which contain toxic or caustic components should only be dried in a fume hood.

- Corrosive: Substances which release corrosive vapors when heated should be tested in small amounts.

- The user assumes responsibility for any damage caused by the use of these types of samples.

EN-2

Application Disclaimer

Moisture determination applications must be optimized and validated by the user according to local regulations. Application specific data provided by OHAUS is for reference purposes only. OHAUS waives all liability for applications based on this data.


2. INSTALLATION

2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.

Select a location that is safe and with adequate ventilation. Fire, corrosive or toxic fumes and other hazards associated with the test samples will require specially prepared locations.

2.2 Installing Components

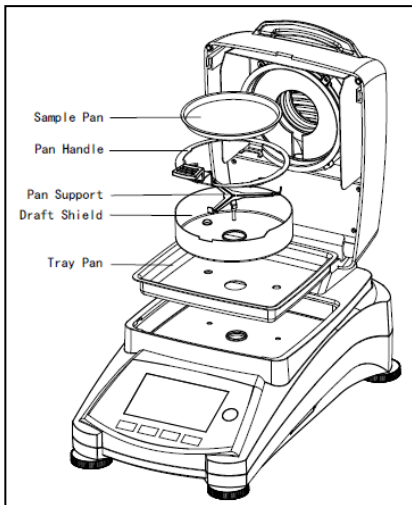2.3 Connecting Power

Connect the AC mains power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

See also section 1, Safety Precautions


After connecting the Analyzer to the AC supply (standby mode), allow the Analyzer to warm up for at least 30 minutes for best results.

Power: On (short press) / **Off / Standby Mode** (long press)

When powered on from standby mode, the Analyzer is ready for immediate use.

2.4 Levelling the Instrument

To level the instrument, adjust the feet so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.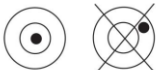


2.5 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

3. OPERATION

3.1 Display and Controls

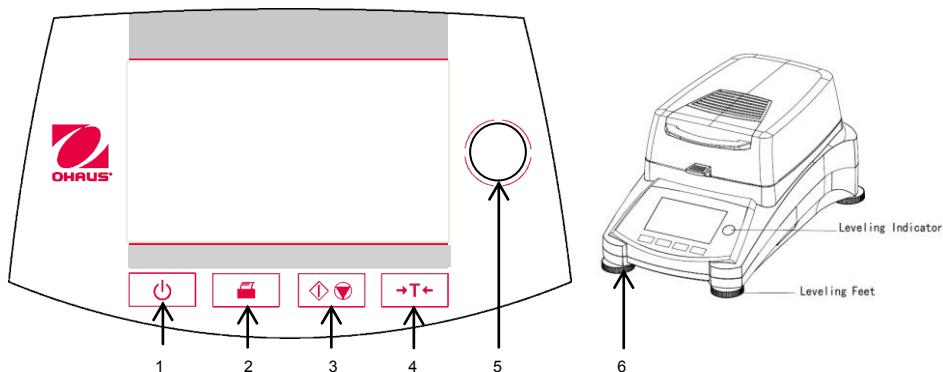
1		Turns On or Off display.
2		When pressed, sends Print command.
3		When pressed, either starts or stops the drying process.
4		When pressed, performs Tare function.
5		Leveling indicator
6	Leveling Feet	Used to level the Moisture Analyzer.

MAKING A SIMPLE MEASUREMENT

If the display is not visible and the unit is plugged into a power source, press the On/Off button.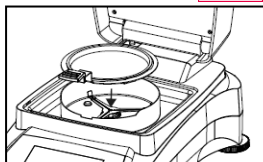


Directly start a measurement.

Open the cover on the Moisture Analyzer. Place the pan handle in the sample chamber. Ensure that the tongue of the pan handle fits exactly in the slot of the draft shield element.

Place the empty sample pan and glass fiber pan in the pan handle.

Note: It is advisable for you to work with the pan handle at all times. The pan handle is ergonomic, safe and provides protection against possible burns due to a hot sample pan.

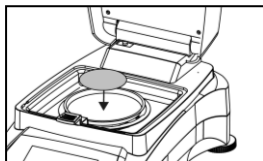Close the cover. This sets the Moisture Analyzer to zero automatically.

Open the cover on the Moisture Analyzer.

Wet the glass fiber pan with a 0.5 to 1 gram of water.

Note: Minimum sample size must be greater than 0.5 gram.

Starting the test

Close the cover, the Moisture Analyzer starts the drying and measurement process automatically.

WARNING: As the pan and sample may still be hot, you should let these cool down before removing the pan from the pan handle!

4. MAINTENANCE

4.1 Cleaning

WARNING: Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

Attention: Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

4.2 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website www.ohaus.com to locate the Ohaus office nearest you.

5. TECHNICAL DATA

The technical data is valid under the following conditions:

Indoor use only

Temperature: 10°C to 40°C

Relative humidity: Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40 °C.

Altitude: Up to 4000 m

Power Input: 100 V – 90 VAC 5A 50/60 Hz or 200 V – 240 VAC 2.5A 50/60 Hz

Voltage fluctuations: Mains supply voltage fluctuations up to ±10% of the nominal voltage

Power load: Max. 450 W during drying process

Power line fuse: 6.3A 250 VAC for 100V-120VAC power supply or 2.5A 250VAC for 200V-240VAC power supply according to the heating element.

Pollution degree: 2

Installation category: Class II

Typical specifications:

Model	MB90	MB120
Capacity (g)	90	120
Readability	0.01%/0.001g	0.01%/0.001g
Temperature Settings	40° C - 200° C	40° C - 230° C
Heat Source	Halogen	Halogen

6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at www.ohaus.com/ce .
	This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .

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P/N 30284480 C © 2016 Ohaus Corporation, all rights reserved.
Ohaus Corporation