
Instruction Manual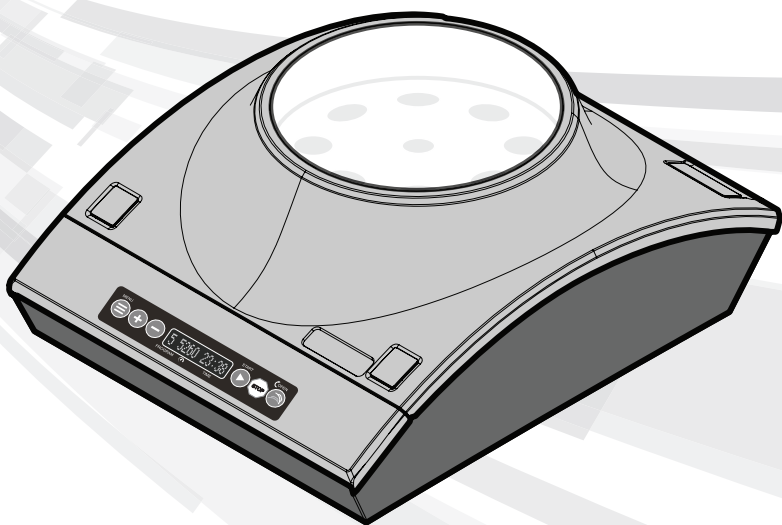

Instruction Manual	2
Manuel d'instructions	21
Manual de Instrucciones	40
Bedienungsanleitung	59
Manuale di Istruzioni	81

Table of Contents

Safety Precautions	3
Technical Data	5
Supplied Equipment	5
Compliance Guidelines	6
EC Declaration of Conformity	7
UKCA Declaration of Conformity	8
Display	9
Rear Panel Switch and Connector	9
Operation	10
Rotor Installation & Removal	11
Centrifuge Rotor Balancing	12
Manual Lid Opening Instructions	12
Centrifuge Instructions	13
Error Messages	15
Troubleshooting	16
Maintenance	17
Warranty	18
Return for Repair	19
Disposal Responsibilities	19

Safety Precautions

This centrifuge is intended for use as a general-purpose laboratory instrument for the separation of components through the application of relative centrifugal force. Applications may include the separation of chemicals, industrial or environmental samples, and bodily fluids such as blood and urine, either alone or after the addition of reagents or other additives. Per Rule 5a of the In Vitro Diagnostics Device Regulations (IVDR 2017/746) this instrument falls within the Class A, non-sterile device designation. This centrifuge must be operated by trained laboratory personnel such as a clinical laboratory technician or technologist or similar in a professional healthcare facility environment.

This instruction manual cannot address all safety hazards. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instrument's suitability to the task.

Caution!

Before starting the device for the first time, please read this instruction manual carefully. This manual contains important information, safeguards, and operating instructions.

1. Always completely inspect the device including the rotor for damage (broken or cracked housing and base, loose fasteners, damage from solvents, etc.) before connecting the power cord and before each use. Do not use this device if any damage is found.
2. Confirm the tubes to be centrifuged fit properly into the rotor, use Heathrow supplied tube adapters and spacers as needed. Do not allow top of tubes to contact inside of lid (scuffing of lid and tube lid damage may occur. Do not allow tubes to fall inside the rotor.
3. Do not attempt to disassemble any part of the centrifuge.
4. Never use this product in any manner inconsistent with these instructions.
5. This product is intended to be used indoors only.
6. Use only the power adapter and grounded (earthed) power cords delivered with the product, use of other power adapters and grounded power cords will void the warranty. Only plug the power cable into properly grounded (earthed) electrical sockets.
7. This centrifuge is intended for separating aqueous solutions in approved test tubes compatible to the RCF of this unit.
8. This centrifuge is not intended to centrifuge very dense materials. Do not use materials with a density greater than 1.2 g/cm³ (1.2 g/mL).
9. Only use rotors supplied by Heathrow. These rotors can withstand 6,000 rpm which is also the maximum speed of the device.
10. Observe general laboratory safety precautions and regulations when using this product.
11. Do not attempt to operate with the cover removed or open.
12. The rotor and lid must always be securely fastened when in operation.
13. Always seat the rotor completely onto the shaft, then hand tighten the rotor nut.

14. Do not use damaged rotors.
15. Do not put hands onto the rotor unless the rotor is completely stopped.
16. Do not move unit while rotor is turning or unit is powered on.
17. Do not immerse the product in water. Injury due to shock or fire may result.
18. The centrifuge rotor must be loaded symmetrically. Operating this unit with an unbalanced rotor will cause severe vibration and damage, which is not covered by warranty.
19. Do not fill tubes while inserted in centrifuging rotors. Spilled liquid may harm unit.
20. Close tube lids before starting centrifuge. Open lids can be torn off during centrifugation and damage the unit. Open lids can cause fluid to be dispersed inside the centrifuge.
21. Do not insert non-standard centrifuge tubes or other foreign articles into the rotor.
22. Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this instrument.
23. Do not use solvent or flammable liquids near this or other electrical equipment.
24. Do not centrifuge or stir flammable, explosive, or corrosive materials.
25. Do not operate in a hazardous or flammable environment.
26. Do not use any solvent on the unit that may attack plastic or cause cracks in the rotors or corrode metal parts in the rotors.
27. Ensure rotors are protected from corrosion and mechanical damage at all times. Rotors must be cleaned with a pH-neutral cleaning liquid and immediately dried.
28. Always work in a manner which endangers neither the user nor any other person.
29. Should the instrument fail to work to specification, immediately stop using it. Clean and troubleshoot the instrument according to the instructions under "Trouble Shooting" before any further use of the instrument. Contact an authorized dealer or the manufacturer if trouble persists.
30. Do not attempt to stop a rotor while unit is running. Doing so may cause the unit to fail and will void the warranty.
31. Only use original manufacturer's rotors, AC adapters, tube adapters and any other spare parts.
32. Repairs are to be performed by trained and authorized service personnel only. Contact Heathrow Scientific® LLC.
33. Opening the instrument housing, rotor or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Heathrow Scientific® LLC for warranty service.
34. Do not attempt to use units that have not been correctly installed or repaired.
35. Do not attempt to disassemble or modify this product.

36. Store at room temperature in a dry area. Do not expose to sunlight, moisture, or extreme temperatures for prolonged periods of time.
37. Caution: Exposure to certain chemicals can damage unit causing corrosion, crazing, and cracking of housing.
38. Only use consumables within their expiration date.
Note: consumables (tubes, reagents, etc.) are NOT INCLUDED with this product.
39. Do not position device in such a way that it is difficult to operate the On/Off power switch.

Warning!

If the unit is not used as recommended by the manufacturer, the overall safety will be impaired.

Warning!

Check MSDS, wear required Personal Protective Equipment, and observe all applicable local and national regulations before dispensing and disposing of hazardous sample.

Warning!

This product does not contain bio-seals as per IEC/EN/CSA 61010-2-020 and cannot provide any level of containment in case of a spill or release of toxic, radioactive, or pathogenic micro-organisms. These materials should not be used with this product.

Warning!

The laboratory procedure should ensure that no person or hazardous substances are present within a 12" (30cm) zone around the device when operating.

Warning!

Products used in vitro in the preparation of samples that have been obtained for examination are considered neither as IVD nor as accessories and fall outside the scope of the Directive unless, based on their characteristics, they are specifically intended for a particular IVD test.

Technical Data

Supplied Equipment (Spare part item numbers):

Note: consumables (tubes, reagents, etc.) to be used within their expiration date and are NOT INCLUDED with this product.

- Centrifuge
- Universal AC power adapter (120779)
- Set of 4 Universal Grounded mains cables for AC adapter for N. America /Japan, UK, EU, and Australia (120875)
- Standard 8-Tube rotor for 15 mL tubes, 6,000 rpm max (120781)
- 1 bag with 8 ea. tube adapters and 24 ea. tube spacers for smaller tubes; rated for up to 6,000 rpm (120783)

Environmental Ratings:

Note: For Indoor use

Altitude: Up to 2,000 m, Atmospheric Pressure: 80kPa – 107kPa

Operating Temperature: 2°C to 40°C

Maximum Relative Humidity:

80% RH up to 31°C decreasing linearly to 50% RH at 40°C

Environmental Storage and Transport Conditions:

Storage temperature: -20°C – 55°C

Atmospheric pressure: 80kPa – 107kPa

Relative humidity: ≤ 80% RH noncondensing

Light exposure to original packaging: no recommendations, machine is enclosed in opaque packaging.

CAUTION: Light exposure on unpacked centrifuge: UV rays reduce the stability of plastics. Do not subject the centrifuge, rotor, or accessories to direct sunlight/UV light.

Precipitation: both unpacked centrifuge and centrifuge in original packaging must be protected from precipitation

Dimensions: 293mm x 366mm x 148mm (11.54”W x 14.41”D x 5.82”H)

Weight: 4.50 kg (9.93 lbs.) (product with empty rotor)

Capacities: Note: Do not use materials with a density greater than 1.2 g/cm³ (1.2 g/mL) 8-Tube rotor: 120 mL total maximum liquid volume

Speed Range:

500 rpm – 6,000 rpm Max. (tolerance: ≤ 2%, typical: +/- 25rpm @ 6,000 rpm), RCF: 33 x g to 4,668 x g, Speed can be adjusted in 1 rpm increments below 1,000 rpm, 5 rpm increments from 1,000 rpm to 6,000 rpm.

Note: Displayed RCF values are for a typical 15 mL conical tube like those offered by Heathrow Scientific LLC. Other tube lengths will result in differing RCF values. Refer to Heathrow Scientific nomograph for approximate RCF values when using other tubes.

Run Time ranges:

Centrifuge mode: 0:30” – 60’:00” or continuous (Note: time increments vary from 0’:01” to 0’:10” to 1’:00 depending on button hold time)

Electrical:

Power Adapter rating: Input: 100 – 240VAC, 2.5A, 50/60Hz

Output: 12VDC, 12.5A, 150W

Input: 12VDC

Current (under load): Approx. 7.7A

Power: Approx. 93W

Warning! ONLY use power cords and power adapter supplied with unit! Do not unplug device while rotor is turning!

Warning! When plugging the power adapter into any outlet/socket. Only use grounded (earthed) electrical outlets/sockets.

Compliance Guidelines:

This mark is the confirmation that the unit conforms to the EU guidelines and has been tested according to the following EU Declaration of Conformity.

Refer to EC Declaration of Conformity for Intended Purpose (IVDR 2017/746)

Conforms to:

IEC/EN 61010-1
IEC/EN 61010-2-020
IEC/EN 61010-2-101
IEC/EN 61326-2-6

ANSI/UL Std. 61010-1
ANSI/UL Std. 61010-2-020
CAN/CSA C22.2 No. 61010-1
CAN/CSA C22.2 No. 61010-2-020

This product has been tested to the requirements of CAN/CSA C22.2 No. 61010-1 3rd edition, including Amendment 1, or a later version of the same standard incorporating the same level of testing requirements.

This product is FDA listed. Product code: JQC; centrifuges for clinical use.

EC Declaration Of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA
European Contact: Emergo, Europe, Princessegracht 20, 2514AP, The Hague, The Netherlands
Product: Clinical Centrifuge
Product Model No(s): 120775, 120776, and variants
Risk Class: Class A, Rule 5a (Article 47, Annex VIII) per IVDR 2017/746

Intended Purpose (IVDR 2017/746)	
What is to be detected or measured?	Not applicable. This centrifuge does not detect or measure any analyte.
Device function such as, screen/ monitor/ diagnose, or aid in diagnosis, prognosis, and prediction:	This centrifuge is intended to separate human samples, including blood, urine and other bodily fluids.
The specific disorder, condition or risk fact of interest that is intended to be detected, defined, or differentiated	Not applicable. This centrifuge is to prepare samples for downstream in vitro diagnostic procedures and it isn't intended to detect or define any clinical condition.
Automation:	Not applicable. This centrifuge is not an automated instrument.
Qualitative, semi-quantitative or quantitative:	Not applicable. This centrifuge does not provide the testing result.
Type of specimen(s):	Human samples, including blood, urine and other bodily fluids.
Testing population:	Not applicable. This centrifuge is for sample preparation and does not provide tests.
Intended user:	This centrifuge should be operated by laboratory professionals only.
For companion diagnostics, the target population and associated medicinal product(s):	Not applicable. This centrifuge is not a companion diagnostics device.

This unit has been constructed and conforms to the following:

Safety Standards and Directives:

EN/IEC 61010-1 ed. 3.1, 2017
EN/IEC 61010-2-020 ed. 3, 2016
EN/IEC 61010-2-101 ed.3, 2018
UL Std. 61010-1 ed. 3, 2019
UL Std. 61010-2-020 ed. 3, 2016
CAN/CSA C22.2 No. 61010-1-12 (R2017) CAN/CSA C22.2 No. 61010-2-020 (R2017)
IVDR 2017/746 (In-Vitro Diagnostic Medical Devices Regulation)
2014/35/EU (Low Voltage Directive)

EMC Standards:

IEC/EN 61326-1: 2020
IEC/EN 61326-2-6: 2020
FCC Part 15, Subpart B: 2015
ICES-003, Issue 7 2012/19/EU (WEEE)
2015/863/EU (RoHS)

Signed:

Gary Kamees
Title: Vice President Product Development/Manufacturing
Location: Vernon Hills, IL
Date of Issue: June 2023

Note: Changes or modifications to the product not expressly approved by the manufacturer could void the user's authority to operate the equipment.

UKCA Declaration Of Conformity

Manufacturer: Heathrow Scientific® LLC, 620 Lakeview Parkway, Vernon Hills, IL 60061 USA

UK Responsible Person: Emergo Consulting (UK) Ltd., Compass House, Viston Park Histon, Cambridge CB24 9BZ

Product: General Purpose Centrifuge

Product Model No(s): 120775, 120776, and and 120xxx (xxx="777-999")

This unit has been constructed and conforms to the following:

Directives and Regulations:

Electrical Equipment (Safety) Regulations 2016

The Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012

The Electromagnetic Compatibility Regulations 2016

EMC Standards and Directives:

EN 61326-1:2020

EN 61326-2-6: 2020

ICES-003, Issue 7

Safety Standards:

EN/IEC 61010-1 ed. 3.1, 2017

EN/IEC 61010-2-020 ed. 3, 2016

EN/IEC 61010-2-101 ed. 3, 2018

UL Std. 61010-1 ed. 3.1, 2017

UL Std. 61010-2-020 ed. 3, 2016

CAN/CSA C22.2 No. 61010-1-12 (R2017)

CAN/CSA C22.2 No. 61010-2-020 (R2017)

Signed:

Gary Kamees

Title: Vice President Product Development/Manufacturing

Location: Vernon Hills, IL

Date of Issue: March 2023

Note: Changes or modifications to the product not expressly approved by the manufacturer could void the user's authority to operate the equipment.


Display

1. MENU Press to step through programming options
2. Plus (+) and Minus (-): to increase or decrease value
3. Start
4. Stop
5. Time Min: Sec or Continuous
6. **OPEN: device will open when rotor is stopped**

Display Status	Possible Cause
COVER OPEN	— Cover is not latched shut.
Program number Speed Time	— Unit is ready to centrifuge.
TILT ERROR	— A sensor in the unit detects device tilt while the rotor is stationary. Only use device on level surface. DO NOT MOVE DEVICE IF POWER IS ON OR ROTOR IS TURNING!!!
SPIN DOWN	— Cycle is slowing down.
ERROR	— An error has occurred. Examine unit for cause. Turn unit OFF then ON to reset program.

Rear Panel Switch & Connector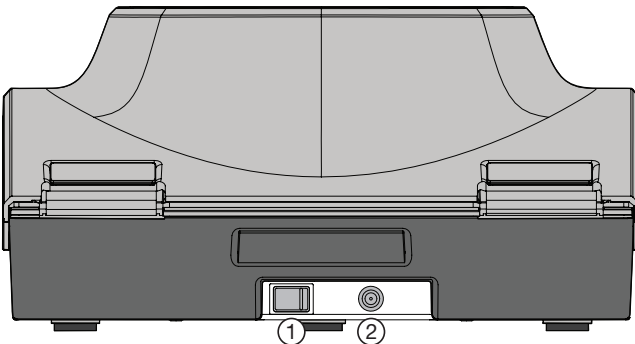

1. Main On/Off Switch
2. Power Input Connector

Operation

1. Remove the contents from the package examining them carefully for breakage, defects or missing parts.
2. Remove the clear protective film from the window on the top of the lid and from the control panel.
3. Place the product on a clean sturdy, level surface at least 12" (30 cm) away from the edge of the counter and other moving equipment. Make sure the AC power adapter and cords are clear of hot surfaces and other hazards. Ensure no person or hazardous substances are within a 12" (30 cm) zone around the product.
4. Make sure the product has ventilation and is not encased in any material that will limit airflow. An overheating situation could occur.
5. Position product so that it is easy to disconnect from mains outlet.
6. Verify that the power switch in the back of the unit is in the OFF position. Plug the DC cord into the rear of the unit and the AC cord an approved outlet.
7. **ONLY** use Heathrow supplied power cord and adapter.
8. **Do not** unplug unit while rotor is in motion.
9. Turn the power switch on the rear of the unit ON. **DO NOT press the start symbol until after removing the cushioning material surrounding the rotor.**
10. Press the open symbol on the control panel to open the lid.
 - Lift the lid and remove the cushion from the top of the rotor.
 - Turn the power switch OFF (set to "O").
 - Unscrew the knurled rotor lock nut by turning it counterclockwise until it disengages from the threaded motor stud.
 - Remove the rotor to access the additional cushioning material around and under the rotor. Discard the cushioning material.
 - See the following ROTOR INSTALLATION & CHANGING section for reinstalling the rotor prior to use.

Prior to every use:

- Inspect unit for cracks, missing pieces, and or any damage that would result in an impaired performance and notify manufacturer BEFORE use.
- Make sure the power switch on the rear of the unit is set to "O" which stands for "OFF".

Rotor Installation & Changing

Warning!

Do not use damaged or cracked rotors!

To remove rotor:

- With one hand, grasp the rotor. Rotors are retained by a central threaded knob. With your other hand, grasp the knob between your fingers and turn the knob counter clockwise ↺ to unscrew the knob until it disengages from the motor shaft.
- Pull up on the center knob firmly. The rotor should easily lift off the motor drive.

—

To install the rotor:

- Check to make sure the two round black O-rings (flexible bands) on the motor are seated in the two grooves of the motor housing.
Fully seated O-rings will slightly protrude from the motor grooves when correctly installed. The O-rings help center the rotor on the motor and keep it from slipping when turning.
- Set the rotor onto the motor housing and push down until seated.
- With one hand, grasp the rotor. Rotors are retained by a central threaded knob. With the other hand, screw the center knob clockwise ↻ until it engages the motor shaft.
- Keep turning the knob until an increase in resistance is felt and it tightens
- **Do not over tighten the knob.**
- **Do not use tools to grasp the knob, damage may result.**


Centrifuge Rotor Balancing

Warning!


Spin balanced loads only! Unbalanced loads may cause damage to the rotor or centrifuge housing from rotor contact.

Do not place the tubes asymmetrically nor load unequal volume tubes. Improper placement of tubes will lead to insufficient centrifugation and may cause serious injury or an accident.


Tubes of equal weight and size should be placed opposite each other. Use additional sample or water in other tubes to provide a balanced rotor. Examples of proper sample balancing are illustrated next.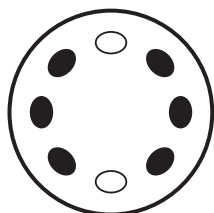

2 Tubes

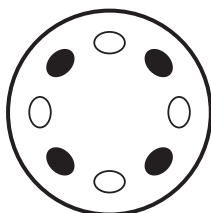4 Tubes

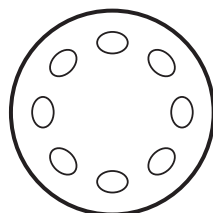
8 Tubes

Warning!

Do not attempt to open lid until rotor has completely stopped! Avoid severe personal injury or property damage from moving parts. Only use rotor compatible with the centrifuge. The lid will unlatch after the cycle has completed.


Manual Lid Opening Instructions

1. Make sure the rotor has come to a full stop and is no longer turning.
2. Turn off power to the unit by using the switch at the rear of the housing.
3. Per international safety standards for centrifuges, the lid can be unlatched by inserting a small tool, such as a screwdriver (not included), into the access hole near the front underside of the centrifuge housing (see arrow). Pushing the tool into the hole (toward the lid) will release the latching mechanism and allow the user to open the lid.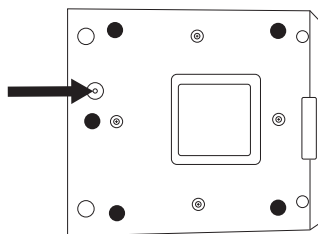

Operating Instructions

1. Only centrifuge tubes containing liquid with a maximum density of 1.2 g/cm^3 (1.2 g/mL).
2. Use ONLY Heathrow Scientific® centrifuge rotors tube spacers and tube adapters when centrifuging. Attempted use of any other rotors will void the warranty and may cause damage to the device or the surroundings. Use of any other rotors may cause INJURY. The Heathrow Scientific centrifuge rotor looks like this:

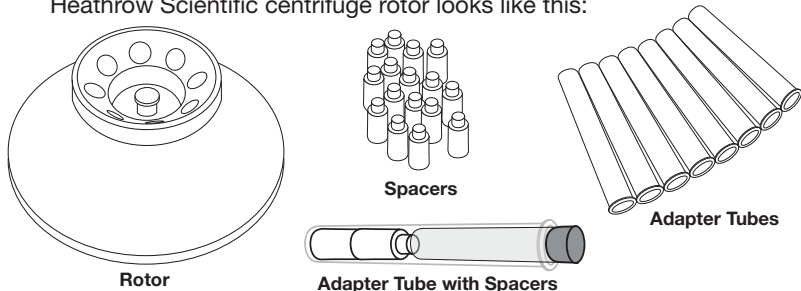3. Make sure power is off on the main power switch on the rear of the housing.
4. Plug the power adapter plug into the connector on the rear of the centrifuge and make sure the blue light on the power adapter is illuminated.
5. Turn on the main power switch
6. The display will immediately light up to display the software/firmware version.

It will then display # SSSSnTTtt where:	
#	Program number from 0 through 9
SSSS	Speed of rotation in rpm or rcf
n	“r” for rpm” or “g” for rcf
TTtt	Time in minutes and seconds or “Cont.” for continuous running.

Note: When first unpacking the centrifuge, press the lid-open symbol to actuate the lid release (you will hear a short clicking sound), open the lid, remove the packing material on top of and under/around the rotor. Install the rotor.

7. Press the lid-open symbol to actuate the lid release (you will hear a click or clicking sound), open the lid and install the rotor. Tighten the retaining knob in the center of the rotor. Close the lid and make sure it's latched. Display will show program number, speed in RPM or RCF, and run time.
8. Press the “+” symbol to select or enter a program number.
Note: Program “0” does not get stored in memory. Programs “1” through “8” can be saved in memory.
Note: Programs “1” through “8” can be overwritten when the rotor is stopped.
Note: When the centrifuge is first turned ON, the last program that was run will be automatically displayed.

9. Press the MENU symbol to select the speed. Either in revolutions/minute (rpm, indicated by the blinking letter “r”) or Relative Centrifugal Force (RCF in “xg” indicated by the blinking letter “g”).
10. Press the MENU symbol again to select speed value. Increase speed value by pressing “+” or “-“ symbols. Press MENU symbol again to save selected value.
11. Time value in “minutes:seconds” or “continuous” (CONT) will now blink. Enter time value and press MENU symbol. Display will now show “+Save Change”. Press “+” to save value or “-“ symbol to change.
12. You have now saved speed, time, and program number.
13. A short press of the START symbol will initiate the program. The rotor will start turning and accelerate to the selected speed, the remaining run time will be displayed. When the remaining run time is 00:00, the rotor will brake to a stop. The centrifuge may slightly overshoot the selected speed momentarily before settling at the programmed value.
14. The time to accelerate to the programmed speed is dependent on the load in the rotor and the selected speed. If the selected time is too brief, the rotor may not achieve the programmed value. When the programmed time is 00:00, the electronic brake slows the unit to a stop. The time to decelerate to a stop is longer than the acceleration time to help prevent the separated fluid components from remixing.
15. When the rotor stops, the centrifuge will emit a “beep” followed by the lid release momentarily activating allowing the lid to open slightly. The phrase “# Complete” will be displayed. Lifting the cover, will change the display to “Cover Open”. Lift the lid upwards and to the rear of the centrifuge to access tubes in the rotor. When completely open, the lid will be tilted slightly rearward and will remain in that position until the operator moves it forward and down.
16. To repeat the existing program, press the lid shut and press the START symbol. Alternately, the user can select a different program by pressing either the “+” or “-“ symbols.

Note: To stop rotor rotation before the timer gets to 00:00 press the STOP symbol. The remaining time will stop counting down and the speed will display the decreasing speed of the rotor.

Warning: do not unplug the unit while the rotor is turning.

Error Messages:

If an error happens, the unit sounds a tone beep and the display will indicate the error.

Error Status	Resolution
Motor Overload	<ul style="list-style-type: none">— Something is interfering with the rotor or the loaded tubes are too heavy. Clear the rotor and reset.
Balance (when centrifuging)	<ul style="list-style-type: none">— Inspect the tubes for equal tube fill or improper placement. Once corrected, rerun.— If the balance error continues to happen, remove the tubes and determine if the balance error still persists with an empty rotor.— Remove/inspect/reinstall rotor.
Temperature	<ul style="list-style-type: none">— The unit has exceeded the normal operating temperature.— Turn off the unit and allow to cool.
Excessive Tilt	<ul style="list-style-type: none">— The unit detects tilt before the rotor starts spinning (do not move unit while rotor is turning)— Make sure the unit is placed on a level surface. Make sure the rotor is undamaged and properly installed. Once corrected, rerun.
Cover Open	<ul style="list-style-type: none">— The lid has unlatched/open. Check for proper operation of the lid latch mechanism. The lid should stay locked during the entire cycle.

Troubleshooting

Trouble	Resolution
No power present	<ul style="list-style-type: none">– Verify that the proper AC input power cord is securely plugged into the power adapter.– Verify that the proper AC input power cord is fully plugged into a powered wall outlet. A blue indicator light will illuminate on the power adapter indicating that it has power.– Verify that the round low voltage output cable of the AC adapter is plugged into the device.– Verify that the power switch on the rear of the device is turned on
Unit or display is not operating normally	<ul style="list-style-type: none">– Turn off the unit, wait 1 minute to allow the internal components to discharge and reset, and then turn power on
Excessive vibration or excessive noise	<ul style="list-style-type: none">– Inspect the tubes for equal fill or improper placement– Inspect the rotor for improper installation– Remove the tubes and determine if the noise persists with an empty rotor– Inspect the housing to lid interface to make sure all lid cushions are present.
Lid will not close and/or latch	<ul style="list-style-type: none">– Verify that nothing is blocking the lid from fully closing– Verify that nothing has fallen into the lock mechanism opening
Lid will not open	<ul style="list-style-type: none">– If for any reason you need to manually open the lid due to an error or power loss, please perform:<ul style="list-style-type: none">– Remove anything placed on top of the device.– Turn off the unit, remove power cord.– Make sure the rotor has stopped completely.– Use a thin rod as a tool (a #1 cross-head type screwdriver works well) and insert it into the opening on the bottom (approx. 0.5" or 13 mm) until you feel resistance.– Press gently but firmly. You will feel a mechanical movement within the unit and the lid latch will release.– Remove the rod, set the unit on the feet– Remove your tubes and reclose the lid

Maintenance

Your centrifuge is normally maintenance-free. Clean unit only when it is not plugged into an electrical outlet. When necessary, the housing and rotor can be wiped using a damp cloth and a mild, non-corrosive detergent.

Warning!

Do not use any solvent on the unit that may attack plastic or cause cracks in the rotor. Ensure the rotor is protected from corrosion and mechanical damage. The rotor must be cleaned with a neutral cleaning liquid

If **hazardous** material is spilled on the unit and or leaked into the unit, do not run the device unless it is safe to do so.

Disassembly of Rotors for Cleaning:

1. Remove rotor from unit
2. Remove screws from the bottom of the rotor
3. Clean as needed
4. Reassemble rotor, making sure that the alignment tab is engaged into the slot.
5. Test run reassembled EMPTY rotor to ensure it is still balanced.
6. **Note:** Ensure all parts are thoroughly dry prior to operation.
7. **Note:** DO NOT autoclave rotor or its parts!

Hazardous Chemicals:

1. Cleaning and decontamination may be necessary as a safeguard before laboratory centrifuges, rotors, and any accessories are maintained, repaired, or transferred. Therefore, follow **ALL** instructions provided in the manual pertaining to this process.
2. In the event hazardous material is spilt, the user is responsible for carrying out the appropriate decontamination procedure.
3. Before using any cleaning or decontamination methods except those recommended by the manufacturer, users should check with the manufacturer that the proposed methods will not damage the equipment.

Warranty

IN NO EVENT WILL HEATHROW SCIENTIFIC® LLC'S OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

Limited Warranty

Heathrow Scientific® LLC warrants that your centrifuge will be free from defects in workmanship and material for five years from the date of purchase.

If you believe that there is a defect in the product, you must, during the warranty period, notify Heathrow Scientific® LLC, provide proof of purchase, and return the product to Heathrow Scientific® LLC with a Return Authorization form. To obtain a Return Authorization form, please call 1-847-816-5070. If Heathrow Scientific® LLC is properly notified and, after inspection, confirms that there is a defect and the warranty period has not expired, Heathrow Scientific® LLC will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, HEATHROW SCIENTIFIC® LLC MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. HEATHROW SCIENTIFIC® LLC WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

- ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (THE CENTRIFUGE IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)
- ANY PRODUCT THAT HAS BEEN, IN HEATHROW SCIENTIFIC® LLC'S SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN HEATHROW SCIENTIFIC® LLC.
- ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.
- ACCESSORY PARTS, SUCH AS RUBBER AND PLASTIC PARTS THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

Return for Repair

Important:

Transporting hazardous materials without a permit is a violation of federal law.

Heathrow Scientific® LLC will not accept any product return that is not appropriately cleaned and decontaminated. In the unlikely event of repair, or when damage to the unit necessitates return, contact Heathrow Scientific® LLC, obtain return authorization, and complete a Heathrow decontamination form BEFORE sending your product for service.

Disposal Responsibilities

The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.

If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact us for more information on the environmental performance of our products.

620 Lakeview Parkway • Vernon Hills, IL 60061
Phone: 1-847-816-5070 • Fax: 1-847-816-5072
Email: info@heatsci.com
Web: www.heathrowscientific.com

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All rights reserved.
©Heathrow Scientific® LLC