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Warranty

In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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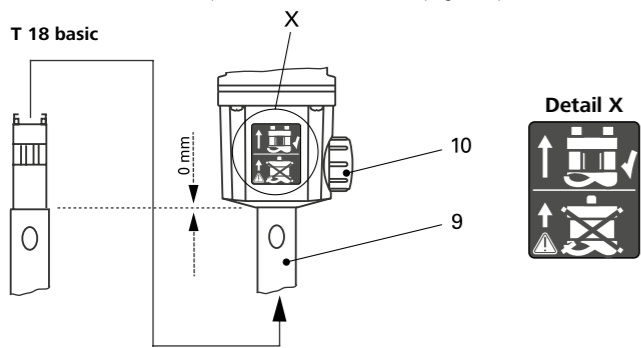
Assembling of the dispersion tool:

- Open the knurled screw (pos. 10) so that the thread does not protrude into the hole.
- Insert the dispersing element (pos. 9) from the bottom into the drive unit as far as the stop. Ensure that the dispersion tool is aligned correctly when inserted (refer to picture: the tabs must project sideways when viewed from the front of the drive unit/dispersion tool). The shaft will audibly click into place when a small amount of force is applied. The dispersion tool is correctly mounted when the edge of the projecting part is aligned with the lower edge of the drive unit (see fig. 2).
- Now secure the dispersing element by screwing the knurled screw tight (pos. 10).

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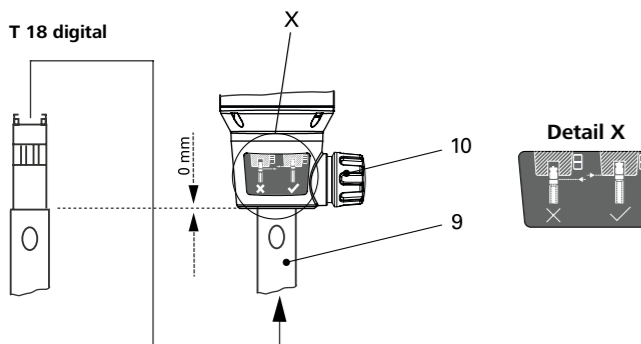
IKA Works Guangzhou
Phone: +86 20 8222 6771
eMail: info@ika.cn

INDIA

IKA India Private Limited
Phone: +91 80 26253 900
eMail: info@ika.in

BRAZIL

IKA Brazil
Phone: +55 19 3772 9600
eMail: sales@ika.net.br

POLAND

IKA Poland Sp. z o.o.
Phone: +48 22 201 99 79
eMail: sales.poland@ika.com

UNITED KINGDOM

IKA England LTD.
Phone: +44 1865 986 162
eMail: sales.England@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com

IKAworldwide

IKAworldwide /// #lookattheblue

@IKAworldwide

Technical specifications may be changed without prior notice.

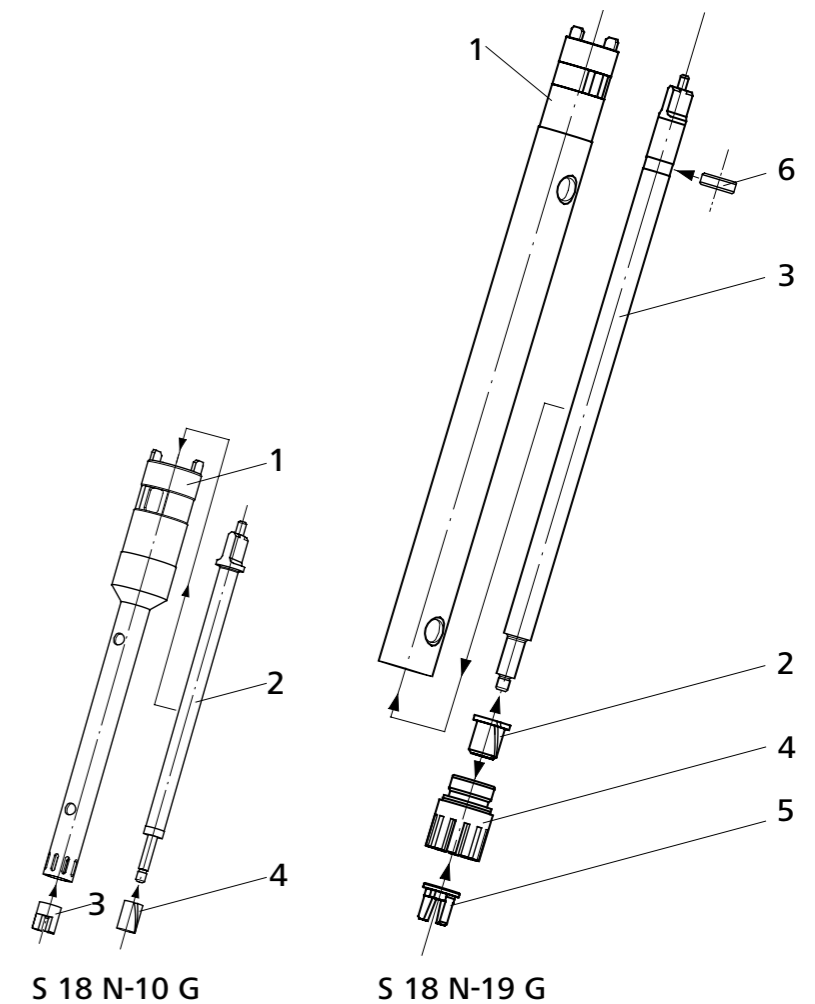
IKA

designed for scientists

Dispergierwerkzeuge/ Dispersion tools/ Outils dispersants

S 18 N - 10 G

S 18 N - 19 G

20000005142c_3567800_S_18_N_082021_print/web

Technische Information
Technical information
Informations Techniche

DE
EN
FR