IKA

designed for scientists

IKA R 2722 H-Stand
IKA R 2723 Telescopic stand

5

/// DEUTSCH

13

/// FRANÇAIS

21

/// ITALIANO

29

/// POLSKI

37

/// TÜRKÇE

45

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9

/// ENGLISH

17

/// ESPAÑOL

25

/// PORTUGUÊS

33

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41

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R 2722

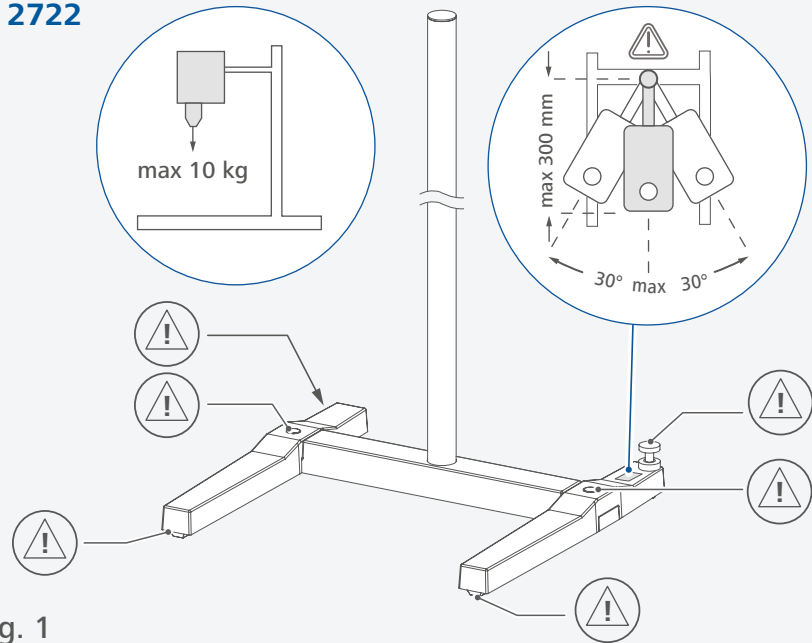
Fig. 1

R 2723

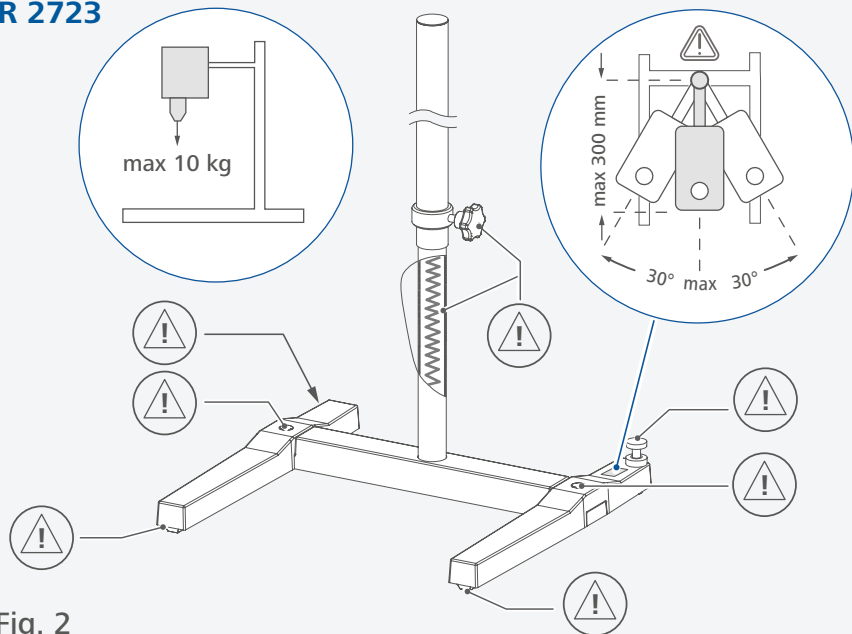
Fig. 2

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	EU Declaration of conformity	9
	Explication of warning symbols	10
	Safety instructions.....	10
	Intended use.....	11
	Unpacking	48
	Assembly	50
	Cleaning	55
	Technical Data	55

Source language: German

EU Declaration of conformity

We declare under our sole responsibility that the product to which this declaration relates is in conformity with directive 2006/42/EC and the following standard or other normative document: EN ISO 12100.

A copy of the complete EU Declaration of conformity can be requested at sales@ika.com.

Explication of warning symbols

/// Warning symbols

Danger! Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.

Warning! Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Caution! Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Notice! Indicates practices which, if not avoided, can result in equipment damage.

/// General Symbols

A — Position number

Correct / Result!
Shows the correct execution or the result of an action step.

Wrong!
Shows the incorrect execution of an action step.

Note!
Displays action steps that require particular attention to detail.

Closed lock!
Indicates the status of "locked".

Open lock!
Indicates the status of "Unlocked".

Safety instructions

/// General information

- › **Read the operating instructions completely before starting up and follow the safety instructions.**
- › Keep the operating instructions in a place where it can be accessed by everyone.
- › Ensure that only trained staff work with the device.
- › Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- › Install the stand on a flat, stable, clean, non-slip, dry and fire-resistant surface.
- › The device must only be used in a technically perfect condition. The correct functioning of the stand must be checked at regular intervals.
- › Prior to each use, always check the stand, stirrer/disperser and accessories for damage. Never use damaged parts.

- › You must also follow the safety instructions contained in the operating instructions for the stirrer or disperser.
- › Check that the stirrer / disperser is installed correctly before use and at regular intervals.

Caution!

- › Check the mounting position of the stirrer / disperser to avoid the danger of the unit tipping over.
- › Observe the maximum permissible load-bearing capacity (see "Technical data").
- › The adjustment screw can be used to compensate for unevenness on the set-up plate. Use the lock nut to lock the adjustment screw.

Notice!

- › Pay attention to the marked sites in Fig. 1 and Fig. 2.

/// Telescopic stand R 2723

Caution!

- › When adjusting the height of the movable stand rod, ensure that nothing becomes caught between the devices or other objects attached to the bracket, e.g. the lower end stop or lip of the container.
- › The stand rod must be held in place when loosening the knurled screw (avoid rapid increases).
- › Always store the stand in a relaxed state.

Intended use

The **R 2722** H-stand and the **R 2723** Telescopic stand are suitable for large, heavy devices and attachments that require absolute stability.

The R 2723 Telescopic stand should only be used to hold, raise and lower stirrers and dispersers. With the pneumatic spring stand rod, instruments and attachments can be raised and lowered smoothly and without difficulty.

The safety of the user cannot be ensured:

- › if the stand is used in conjunctions with accessories not made or recommended by the manufacturer
- › if the stand is not used for the intended purpose as specified by the manufacturer.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited