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Warranty

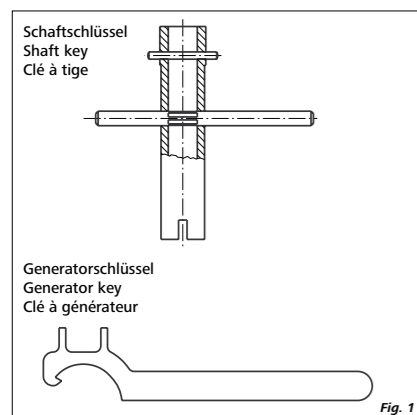
In accordance with IKA® warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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
The tool kit is part of the delivery scope of the disperser.

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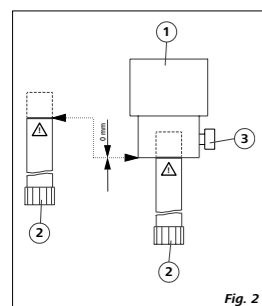Assembling the dispersion tool (overview Fig. 2)

- Open the knurled screw (item 3) so that the thread does not protrude into the hole.
- Insert the dispersing element (item 2) into the drive unit (item 1) as far as the stop. After some resistance (depending on the model of the drive - pressure ball) you will hear the shaft clicking into place. The dispersing element is fitted correctly if the recessed edge of the dispersing element is flush with the bottom edge of the drive unit.

- **Attention:** Now secure the dispersing element by screwing the knurled screw tight (item 3).

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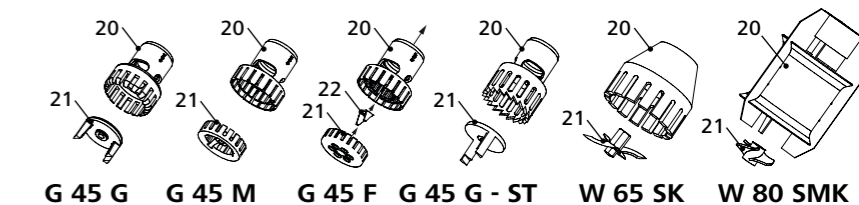
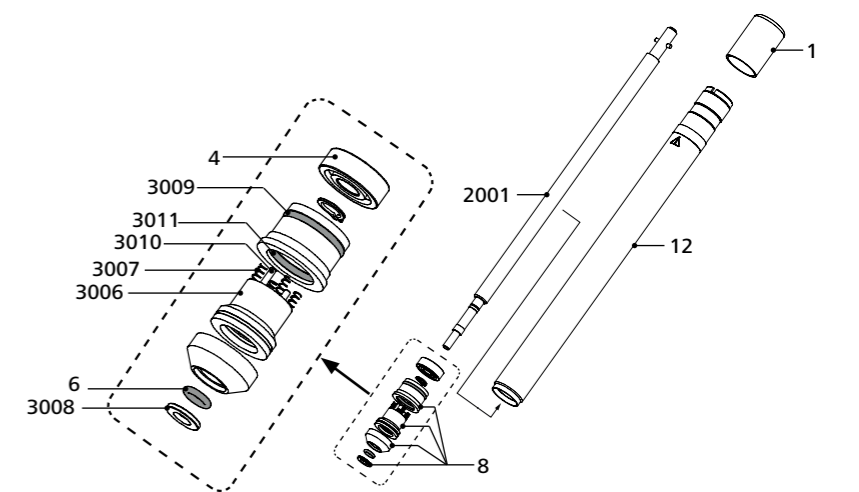
S 50 KG - HH - W 65 SK

S 50 KG - HH - W 80 SMK

Discover and order the fascinating products of IKA online:

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<p> Safety instructions</p>
<div><div></div><div></div></div> <ul style="list-style-type: none">The dispersion tool may heats up during operation. In particular in the processing of hot medium, the tool must be cooled before it can be removed from the drive.When in operation the dispersion tool can become hot. If the dispersion tool is not inserted into the drive flange correctly or the medium is hot, it can become extremely hot. <div> <ul style="list-style-type: none">Don’t touch rotating parts during operation. Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. There is a risk of splashing and evaporation of liquids; body parts, hair, clothing and jewelry getting caught. The work has to be interrupted immediately if you notice unusual noise and/ or increasing emission of liquid from the upper, side-wise rinsing drilling. The bearings must be controlled, and replaced if necessary. Note the operating instructions of the disperser unit. Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on. Do not use any damaged dispersion tools. <i>Attention!</i> Never run the dispersion tool dry. Without cooling by the medium the sealing will be destroyed. The dispersion tool is not suitable for continuous operation. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. Usually a few seconds are sufficient, maximum duration is one minute. Longer application times bring no improvements, increase however the sample temperature substantially. Dispersing causes the medium to heat up. Note the min./max. immersion depth in the “Technical data”. The function of the dispersion tools depends on the condition of the sharp edges of rotor and stator (integrated in the shaft tube). Abrasive media can round off these edges fast, whereby the disperse effect diminishes. </div>
<p>Product information</p>
<div>The bearing of the S 50 KG-HH - shaft consists of a ball bearing and a oneway friction ring seal. The friction rings are made of carbide. The generator (dispersion head) consists of a rotor and a stator. These parts can be unscrewed from the shaft. Note that rotors and stators form matching pairs and must not be confused.</div>
<p>Application instructions</p>
<div>Application: Wet crushing, suspension, emulsifying (Batch-operation).</div>
<p>Spare parts list</p>
<div> <div>Spare parts diagram see front page.</div> <div>For orders of spare parts, please specify the type of dispersion tool and the designation of the spare part.</div> </div>

Item	Designation	Item	Designation
1	Protection cap	3007	Sping
4	Grooved ball bearing	3008	Washer
6	O-ring	3009	O-ring
12	Shaft tube	3010	Pin
2001	Rotary shaft	3011	O-ring
3006	Axial face seal		

Maintenance and cleaning

Cleaning:

For pre-cleaning the dispersion tool, operate it in a detergent for dissolving the residuary of the substance. Due to the large flow rate, rotor and stator are cleaned to a large extent. The dispersion tool must be divided and cleaned immediately after working, so that the adhering residuary of the substance don’t cultivate unwanted bacterial cultures.

For cleaning purposes disassemble the dispersion tool as described below by the shaft and the generator key, both included in the tool kit of the dispersion drive:

- Unscrew the rotor **(21)** and blade **(22)** from the rotary shaft **(2001)** using the generator and shaft wrenches included in the tool kit. The generator wrench is used to hold the rotor while unscrewing it by turning the shaft from above with the shaft wrench. ***Caution:*** The dispersion head **G 45 G** and **G 45 G - ST** has two prominent blades.
- Unscrew the stator **(20)** from the shaft tube **(12)** (left-hand thread). Insert the cross pin (3 mm) of the shaft wrench into the slot on the end face of the shaft tube. The stator can now be unscrewed by turning it in the clockwise direction with the generator wrench.
- In this position the components of the slide ring sealing **(8)** can be pulled off the rotary shaft **(2001)**.
- Press the rotary shaft **(2001)** with ball bearing **(4)** down and out of the shaft tube.
- The slide ring sealing **(8)** must be replaced if they are worn. The friction surfaces of the sealing must be clean.

Reassemble the dispersion tool in the reverse sequence.

Note: when assembling the slide ring sealing **(8)** to the shaft, first assemble the sealing without the o-ring **(6)**, then assemble the o-ring **(6)** until to the groove, then assemble the washer **(3008)** and the re rotor **(21)**.

Attention! The rotor must not be tightened too strongly, otherwise the thread will be damaged.

Repair:

Please send the dispersing tool for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For repair, please request the **“Decontamination Certificate”** form **IKA**® or use the download printout of it from **IKA**® website: **www.ika.com**. If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

<p>Technical Data</p>							
		S50KG-HH-G45G	S50KG-HH-G45M	S50KG-HH-G45F	S50KG-HH-G45G-ST	S50KG-HH-W65SK	S50KG-HH-W80MSK
Working range	l	0.5–20	0.5–15	0.25–10	0.5–20	1–10	1–50
Stator/Rotor (Ø)	mm	45 / 36	45 / 40.5	45 / 40	45 / 38	65 / 55	80 / 34
Gap between rotor and stator	mm	0.5	0.25	0.5	0.5	-	-
Max. allowable speed	rpm	10000					
Max. circumferential speed	m/s	18.8	21.2	20.9	19.9	28.8	17.8
Min./max. immersion depth	mm	70 / 260	70 / 250	70 / 250	70/260	80–350	140–350
Materials in contact with medium		FFPM, WC, AISI 316 L					
pH range		2–13					
Suitable for solvents		Yes					
Suitable for abrasive substances		No					
Max. working temperature	°C	150					
Sterilization methods		wet chemical					
Working range pressure	bar	6					
Working range vacuum	mbar	1					
Ultimate fineness, suspensions	µm	40–100	25–50	10–30	40–100	50–200	-
Ultimate fineness, emulsions	µm	10–30	5–20	1–10	10–30	5–15	-

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