

# Amendment

## Amendment to XDS5/10 and 35i Silencer Kit (A505-97-880 Issue C)

The following section has been updated:

### 3 INSTALLATION

#### 3.2 Install the XDS silencer

Ensure the pump is switched off before fitting.

1. Clean the face of the pump exhaust with a lint free cloth.
2. Place the centring ring (Figure 1, item 5) in the end of the NW25 flange and seat the silencer in the pump.

**Below is an additional comment that needs to be added to the instruction manual**

*Note: Exhaust hole should be pointed downward and away from the operator to avoid any foreign body, process material or condensate that may be evacuated on start up.*



This page has been intentionally left blank.