IKA

designed for scientists

IKA Rotastand

5
/// DEUTSCH

13
/// FRANÇAIS

21
/// ITALIANO

29
/// PORTUGUÊS

37
/// TÜRKÇE

9
/// ENGLISH

17
/// ESPAÑOL

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33
/// POLISH

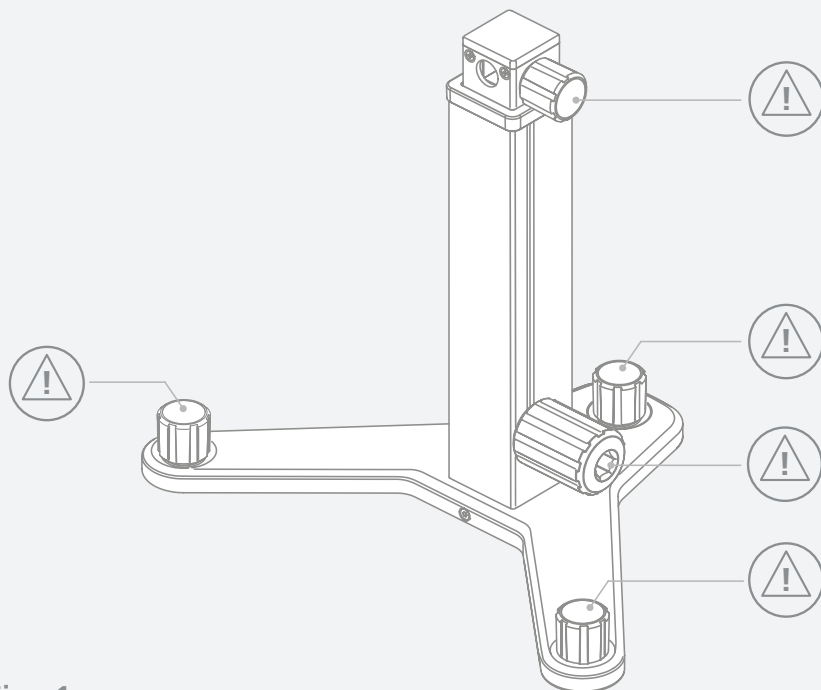
Fig. 1

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	EU Declaration of conformity	9
	Explication of warning symbols	9
	/// Warning symbols	9
	/// General Symbols	10
	Safety instructions.....	10
	/// General information.....	10
	/// Device setup.....	10
	/// Working with the device.....	10
	/// Accessories.....	10
	Intended use.....	10
	/// Use.....	10
	/// Area of use (only indoors).....	11
	Unpacking	11
	/// Unpacking	11
	/// Scope of delivery.....	11
	Assembly	11
	Maintenance and cleaning	12
	Technical Data	12
	Warranty	12

Source language: German

EU Declaration of conformity

We declare under our sole responsibility that this product corresponds to the directive 2006/42/EC and conforms with the following standard or normative document: EN ISO 12100.

A copy of the complete EU Declaration of conformity can be requested at sales@ika.com.

Explication of warning symbols

/// Warning symbols

DANGER

Indicates an (extremely) hazardous situation, which, if not avoided, will result in death, serious injury.

WARNING

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

CAUTION

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

NOTICE

Indicates practices which, if not avoided, can result in equipment damage.

/// General Symbols

A — Position number
Indicates device components relevant to actions.

Safety instructions

/// General information

- › **Read the operating instructions completely before starting up and follow the safety instructions.**
- › Keep the operating instructions in a place where it can be accessed by everyone.
- › Follow the safety instructions, guidelines, occupational health, safety and accident prevention regulations.

/// Device setup

- › Set up the device in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- › All screw connections must be properly tightened.
- › Check the mounting position of the viscometer to avoid the danger of the unit tipping over.

/// Working with the device

Notice!

- › Risk of tipping exists when tilted the stand at an angle of being over 10° with full load.
- › Observe the maximum permissible load-bearing capacity (see "Technical data").
- › When moving the stand, ensure the device is at the lowest end stop position.
- › When used as intended and in accordance with the operating instructions, there is no danger to people, the environment or the device. The manufacturer cannot be held responsible for any damages as a result of incorrect use of the device or non-compliance with the safety instructions.

/// Accessories

- › Protect the device and accessories from bumping and impacting.
- › Check the device and accessories beforehand for damage each time when you use them. Do not use damaged components.

Intended use

/// Use

- › The IKA Rotastand has been designed for holding, lifting and lowering IKA Rotavisc.

/// Area of use

- › Indoor environments similar to that a laboratory of research, teaching, trade or industry area.
- › The safety of the user cannot be guaranteed:
 - If the device is operated with accessories that are not supplied or recommended by the manufacturer.
 - If the device is operated improperly or contrary to the manufacturer's specifications.
 - If the device is modified by the third parties.


Unpacking

/// Unpacking



- › Unpack the device carefully. Any damage should immediately be reported to the carrier (mail, rail or freight forwarding company).

/// Scope of delivery

M	Rotastand pillar
N	Rotastand base
O	Screw washer
P	Hexagon socket screw
Q	Hexagon socket screw
R	Screwdriver (AF = 2.5 mm) DIN ISO 2936
S	Screwdriver (AF = 5 mm) DIN ISO 2936

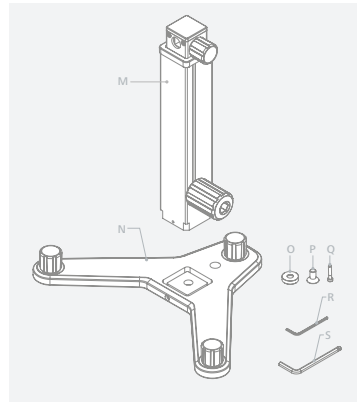Assembly

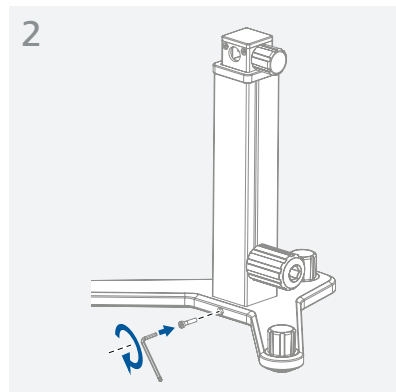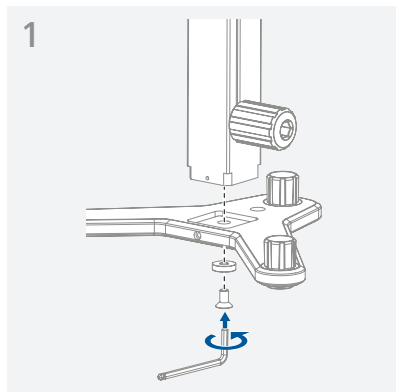Maintenance and cleaning

/// Proper handling

- › The device is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Cleaning:

- › Use only cleaning agents which have been approved by IKA to clean IKA devices. These are water (with tenside) and isopropanol.
- › Wear protective gloves during cleaning the devices.
- › Do not allow moisture to get into the device when cleaning.
- › Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Ordering spare parts:

- › When ordering spare parts, please give:
 - Device type,
 - Position number and description of spare part, see www.ika.com.

Repairs:

- › Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.
- › For this, use the **"Safety Declaration (Decontamination Certificate)"** form which you can obtain from IKA or can download a version for printing from the IKA website at www.ika.com.
- › If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

Technical data

Stroke max.	mm	200
Diameter (Ø)	mm	16
Dynamic load	kg	5
Dimensions (W x D x H)	mm	351 x 246 x 318
Weight	kg	4.8

Warranty

- › In accordance with IKA warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.
- › The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited