

**Tuttnauer** Lab

Sterilization & Life Science Solutions

# T-Lab Eco Line

Affordable Performance and Trusted Reliability  
for your laboratory sterilization needs

LEADING IN  
STERILIZATION  
SINCE  
**1925**

Laboratory



# The Economic Laboratory Autoclave From The World's Premium Manufacturer

Vertical Autoclave



Economic Line



Reliable Performance



Minimal Laboratory Resource Consumption



**T-Lab Eco** was developed especially for educational institutions and research facilities. Designed with simplicity and economy in mind.

This vertical floor-standing autoclave covers the fundamental needs for general lab sterilization with the aim of increasing the **productivity** of your laboratory.

Offering the best universal capacity, together with the **optimization** of resources such as water, power and operating time results.

An affordable and reliable solution you can count on for managing your laboratory **workload**.

## Sterilization Application

Intended for the sterilization of a wide range of liquids and solids such as glassware, metal utensils and other laboratory items.

This reliable workhorse is ideal for the general laboratory in need of easy operation, safety and cycle features for day-to-day needs.

### LABORATORY APPLICATIONS



### SOLID, GLASSWARE LOADS



# Making Your Process Easier, Accurate and Affordable

Intuitive, user friendly technology, including colour touchscreen and multiple programmes for sterilization.

Feature Advantages Include:

## Reliable Performance

**This basic product will do what you need and expect**

- Solid and glassware loads
- Liquid loads
- Efficient sterilization time including cooling
- Two PT100 sensors as standard

## Affordable

**Less in More**

- Built to last without paying for the features and parts you do not need

**Greater Value at Lower Cost**

Less water wasted, less electricity consumed, less time spent on maintenance and repair.

## Plug & Play

**This autoclave's plug and play mode does not require dedicated installation connections, simply a power source**

- Multicolour user-friendly touch screen
- Connectivity to R.P.C.R
- Stores the last 1000 cycles in built-in memory
- Portable - comes standard with lockable wheels

## Versatile

**Adaptable for All Your Basic Needs**

- Sterilizes liquids, solids, sealed and unsealed bottles
- Performance and quality within a small space
- Cylindrical autoclaves – easier to operate, less expensive to maintain, and more reliable over a longer working life

## Safety

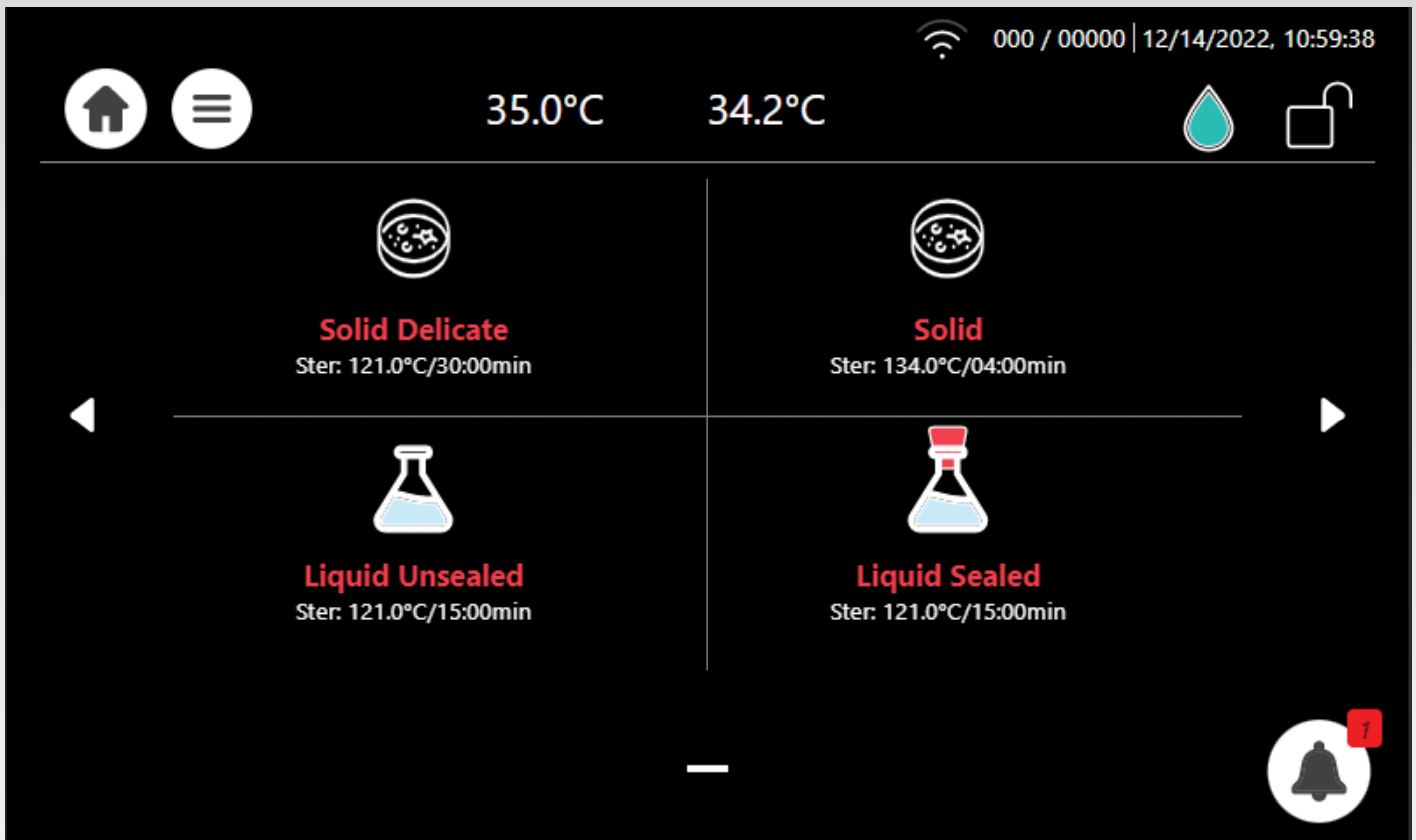
**Meets International Regulations**

- International standards + PED certified





EQUIPPED WITH UNPARALLELED FEATURES, UNIQUE TO T-LAB ECO

- Two PT100 sensors
- Built-in cooling fan
- High-resolution touchscreen
- Built-in memory: 1000 cycles





## STERILIZATION PROGRAMMES

Cycle icon	Cycle name	Temperature	Sterilization time (minutes) V60	Sterilization time (minutes) V85	Load Type	Type of Use
	Liquid Sealed	121°C	15	15	Container with Liquid	For storage
	Liquid Unsealed	121°C	15	15	Container with Liquid	Immediate use only
	Solid	134°C	4	4	Solid Load	Immediate use only
	Solid Delicate	121°C	30	30	Solid Load	Immediate use only

# LOADING CAPACITIES




ERLENMEYER FLASKS Size	V60 Total Units : #Unit x #Basket	V85 Total Units: #Unit x #Basket	BOTTLES / Schott Duran Size	V60 Total Units : #Unit x #Basket	V85 Total Units: #Unit x #Basket
<b>250 ml</b> Ø 85 x 143 mm	13 x 2	13 x 2	<b>250 ml</b> Ø 70 x 143 mm	19 x 2	19 x 2
<b>500 ml</b> Ø 105 x 183 mm	8 x 2	8 x 2	<b>500 ml</b> Ø 86 x 182 mm	12 x 2	12 x 2
<b>1000 ml</b> Ø 131 x 230 mm	5 x 2	5 x 2	<b>1000 ml</b> Ø 101 x 230 mm	8 x 2	8 x 2
<b>2000 ml</b> Ø 166 x 280 mm	3 x 1	3 x 1	<b>2000 ml</b> Ø 136 x 260 mm	4 x 1	4 x 1

\* Loading capacities in this tables are non-binding guidance.

\* Total units per basket have been calculated using International Standard Erlenmeyer's Flasks and Schott Duran.



## Stainless Steel Wire Baskets

2 Stainless Steel wire baskets included	Stainless steel basket - Optional	Leakage tray - Optional
		

## Baskets dimensions

Specifications	V60	V85
Basket Ø x H (mm)	362 x 190	362 x 280
Pail Ø x H (mm)	350 x 320	350 x 450
Tray Ø x H (mm)	345 x 40	345 x 40

## Optional Thermal Printer

### T-LAB Printer – Desktop Direct Thermal Bar Code Printer

- High quality double-walled clamshell design
- 127 mm (5") OD media capacity
- Up to 127 mm (5") per second print speed
- Easy media loading
- Head open sensor
- RS232 connection



Technical Features	
Resolution	8 dots/mm (203 DPI)
Max. print speed	127 mm (5")/second
Max. print width	54 mm (2.13")
Max. print length	2,286 mm (90")
Overall Dimensions	W x H x D (mm) W x H x D (Inch)
	109 x 171 x 209 4.29" x 6.73" x 8.23"
Weight	1.2 kg (2.65 lbs)
Label roll capacity	127 mm (5") OD
Memory	4 MB Flash memory   8 MB SDRAM   microSD card for flash memory expansion, up to 4 GB
Interface	RS-232   USB 2.0   Bluetooth (user option)
Accessories	Windows labelling software CD disk   External universal switching power supply

## Regulations

### **Company Standards and Legislation**

EMC Directive 2014/30/EU

ISO 9001 Quality Management System

2006/42/EC Machinery Directive

2014/35/EU, Low Voltage Directive (LVD)

2012/19/EU WEEE and EU 2017/2102 Restriction of the use of certain hazardous substances (RoHS Directive)

### **Product Standards**

Pressure Vessel – PED 2014/68/EU

IEC 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use –

Part 1: General requirements

IEC 61010-2-040: Safety requirements for electrical equipment for measurement, control, and laboratory use –

Part 2-040: Particular requirements for sterilizers and washer-disinfectors used to treat medical materials

EN 61326-1: Electrical equipment for measurement, control, and laboratory use – EMC requirements –

Part 1: General requirements



# TECHNICAL SPECIFICATIONS

Specifications	V60	V85
Chamber Volume (Litres)	60	85
Weight (kg)	111	120
Chamber Dimensions Ø x D (mm)	390 x 505	390 x 700
External Dimensions W x H x D (mm)	645 x 985 x 810	645 x 985 x 810
Shipping Dimensions W x H x D (mm)	820 x 1250 x 900	820 x 1250 x 900
Ratings	Voltage: 220-240 VAC Single Phase Power: 3000 W Amperage: 13.1 A	Voltage: 220-240 VAC Single Phase Power: 5400 W Amperage: 22.3 A
Wastewater tank	3 Litres	3 Litres
Cooling water tank	6 Litres	6 Litres
Maximum Allowable Working Pressure	2.8 bar	2.8 bar



Your lab infection control success is backed more than a century of global leadership experience

A leader in infection prevention

Products distributed to clinics, hospitals, and research institutions in more than

since  
**1925**

**140**  
countries

Laboratory

**International Sales and Marketing**  
E-mail: [info@tuttnauer.com](mailto:info@tuttnauer.com)  
[www.tuttnauer.com](http://www.tuttnauer.com)

**Tuttnauer Europe B.V.**  
Hoeksteen 11, 4815 PR  
PO Box 7191, 4800 GD Breda  
The Netherlands  
Tel: +31 765 423 510  
Fax: +31 765 423 540  
E-mail: [info@tuttnauer.nl](mailto:info@tuttnauer.nl)

**Tuttnauer USA Co.**  
25 Power Drive,  
Hauppauge, NY 11788  
Tel: +800 624 5836, +631 737 4850  
Fax: +631 737 0720  
E-mail: [info@tuttnauerUSA.com](mailto:info@tuttnauerUSA.com)

Distributed by:



Learn from our experts  
Join our blog [tuttnauer.com/blog](http://tuttnauer.com/blog)