20000005489

**IKA®**

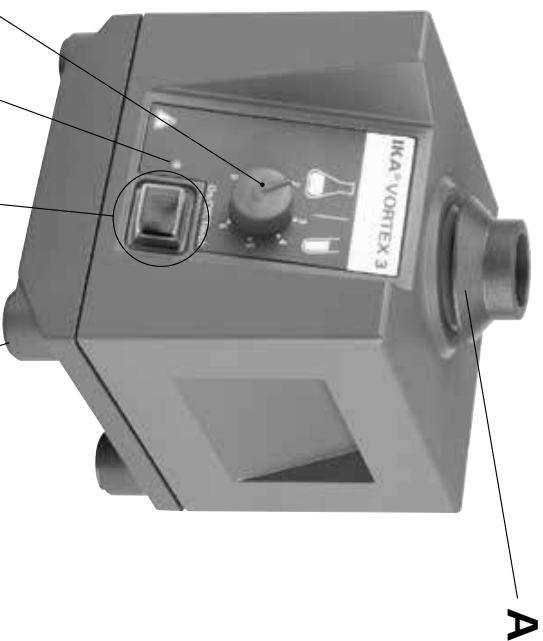
# IKA® Vortex 3

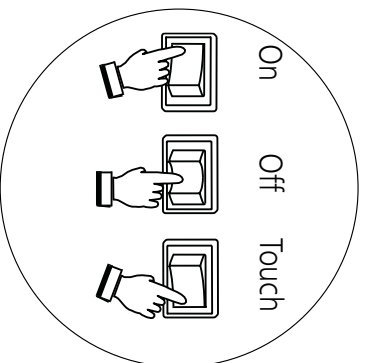
Betriebsanleitung	DE	3
Ursprungssprache		
Operating instructions	EN	9
Mode d'emploi	FR	15
Indicaciones de seguridad	ES	21
Veiligheidsinstructies	NL	22
Norme di sicurezza	IT	23
Säkerhetsanvisningar	SV	24
Sikkerhedsanvisninger	DA	25
Sikkerhetsanvisninger	NO	26
Turvallisuusohjeet	FI	27
Instruções de serviço	PT	28
Wskazówki bezpieczeństwa	PL	29
Bezpečnostní upozornění	CS	30
Biztonsági utasítások	HU	31
Varnostna navodila	SL	32
Bezpečnostné pokyny	SK	33
Ohutussjuhised	ET	34
Drošības norādes	LV	35
Saugos reikalavimai	LT	36

IKA-Motite, Germany  
Reg. No. 004343

**D** Schalter/ Switch/ Commutateur

**Fig. 1**

Pos.	Bezeichnung	Item	Designation	Pos.	Désignation
A	Standardaufsatz V.3.1	A	Standard attachment V.3.1	A	Support standard V.3.1
B	Drehknopf	B	Rotating knob	B	Bouton rotatif
C	Kontrollleuchte	C	Control light	C	Lumière contrôle
D	Geräteschalter	D	Switch	D	Commutateur
E	Gerätefuß	E	Base	E	Pied

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**DE**

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	Page
Declaration of conformity	2
Explication of warning symbols	9
Safety instructions	9
Unpack	10
Correct use	10
Commissioning	11
Accessories	12
Maintenance	13
Warranty	13
Technical data	14

### Explication of warning symbols

Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates practices which, if not avoided, can result in equipment damage.

- For your protection:
- **Read the operating instructions in full before starting up and follow the safety instructions.**
  - Keep the operating instructions in a place where they can be accessed by everyone.
  - Ensure that only trained staff work with the appliance.
  - Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
  - Socket must be earthed (protective ground contact).

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of:

- splashing liquids
- projectile parts
- body parts, hair, clothing and jewellery getting caught.

Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:

- Keep the sample container as vertical as possible.
- Only exert the minimum pressure necessary for the mixing process.
- Use attachments for multiple samples when there is a large number of samples.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- Position the knob (B) at the left stop before starting up. Gradually increase the speed.
- If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.

Reduce the speed if:

- the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the appliance begins to move around because of dynamic forces.
- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.
  - Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
  - Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
  - The appliance is not suitable for manual operation.
  - Sharp edged vessels cause wear debris on the attachments.

Beware of the risk of:

- flammable materials,
  - glass breakage as a result of mechanical shaking power.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
  - **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
  - Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.
  - Always disconnect the plug before fitting accessories.
  - The appliance starts up again automatically following a cut in the power supply.
  - The appliance may heat up when in use.

For protection of the equipment

- The voltage stated on the nameplate must correspond to the mains voltage.
- Protect the appliance and accessories from bumps and impacts.
- The appliance may only be opened by experts.

## Unpack

- **Unpack**
  - Please unpack the device carefully
  - In the case of any damage a fact report must be sent immediately (post, rail or forwarder)
- **Delivery scope**
  - Shaker **Vortex 3**
  - Standard attachment V 3.1
  - Operating instructions

## Correct use

- **Use**
    - ▶ For mixing liquids
    - Touch mode for single test tubes
    - Continuous mode for single or several vessels
  - **Range of use (indoor use only)**
    - Laboratories
    - Pharmacies
    - Schools
- This device is suitable for use in all areas except:
- Residential areas
  - Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed if the appliance is operated with accessories that are not supplied or recommended by the manufacturer or if the appliance is operated improperly contrary to the manufacturer's specifications.

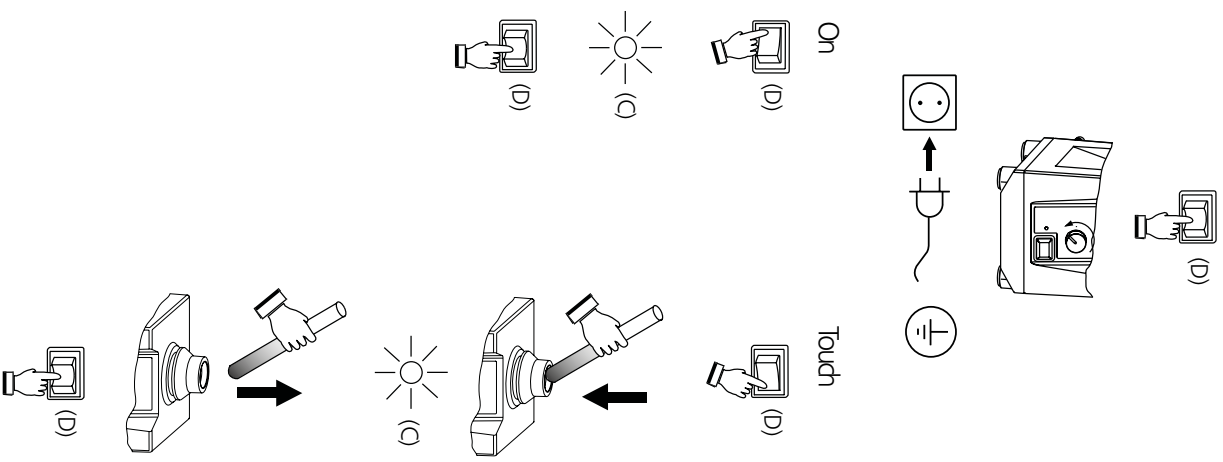
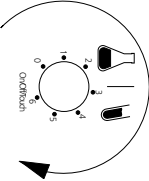
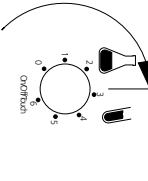
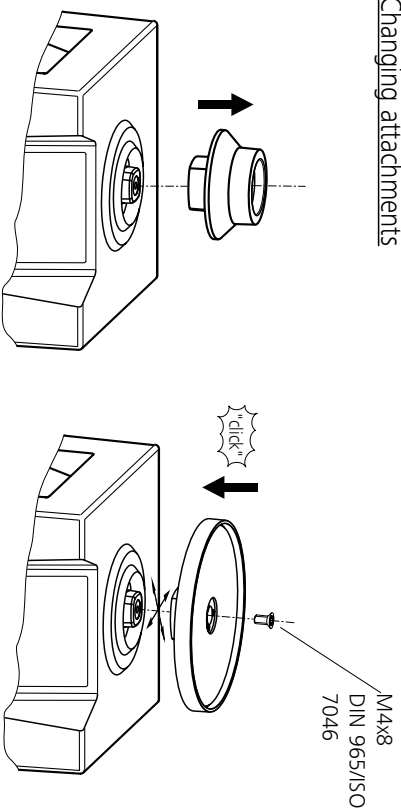
Fig. 2

## Accessories

Possible uses and permitted speed ranges of attachments:

Designation	Description	Id. - No.	Touch Mode	Continuous Mode	Range of speed
 V 3.1 Standard attachment	For test tubes and small vessels upto diameter 30 mm	33 412 00	X	X	
 V 3.2 One-hand attachment	88 mm-plate attachment with rubber pad	33 423 00	X	X	
 V 3.3 Universal attachment	150 mm-plate attachment with rubber pad	33 424 00	-	X	
 V 3.31 Test tube insert	For 54 Eppendorf tubes for V 3.3	33 443 00	-	X	
 V 3.32 Test tube insert	For 18 test tubes 10 mm, for V 3.3	33 439 00	-	X	
 V 3.33 Test tube insert	For 12 test tubes 12 mm, for V 3.3	33 440 00	-	X	
 V 3.34 Test tube insert	For 8 test tubes 16 mm, for V 3.3	33 441 00	-	X	
 V 3.35 Test tube insert	For 8 test tubes 20 mm, for V 3.3	33 442 00	-	X	
 V 3.36 Erlenmeyer flask attachment	For 1 Erlenmeyer-/ round-bottomed flask 100-250 ml, for V 3.3	33 421 00	-	X	
 V 3.37 Microtiter plate insert	For 1 microtiter plate, für V 3.3	33 444 00	-	X	

## Changing attachments

## Using inserts

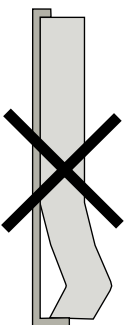
Fig. 4

## Maintenance

The appliance is maintenance-free.

### Cleaning

 →  →  Only use cleansing agents which have been recommended by **IKA®**

Use to remove:

- |                        |   |
|------------------------|---|
| Dyes                   | isopropyl alcohol                           |
| Construction materials | water containing tenside/ isopropyl alcohol |
| Cosmetics              | water containing tenside/ isopropyl alcohol |
| Foodstuffs             | water containing tenside                    |
| Fuels                  | water containing tenside                    |
- Do not allow moisture to get into the appliance when cleaning
- Wear protective gloves during cleaning the devices.

- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

- The rubber of the attachments' pads (V 3.2 and V 3.3) is a natural product, so talcum can discharge. Clean the rubber pad with glycerin.

### Spare parts order

When ordering spare parts, please give:

- Machine type
- Manufacturing number, see type plate
- Item and designation of the spare part, see [www.ika.com](http://www.ika.com), spare parts diagram and spare parts list

### Repair

**Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.**

For this, use the "certificate of compliance" form which you can obtain from **IKA®** or can download a version for printing from the **IKA®** website at [www.ika.com](http://www.ika.com).

If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

## Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## Technical data

Operating voltage	<b>VAC</b>	100 – 120 ± 10%
Design frequency	<b>Hz</b>	60
Nominal voltage	<b>V/Hz</b>	115/ 60
Input power	<b>W</b>	60
Output power	<b>W</b>	10
Speed	<b>1/min</b>	0/ 500 – 2500 infinitely adjustable
Speed adjustment		analog (infinitely) rotating knob in front of device Scale 0 – 6 Shaded pole motor
Speed display		
Drive		4
Agitation stroke	<b>mm</b>	horizontal
Shaking motion		approximately-circular
Perm duration of operation	<b>%</b>	100
Perm ambient temperature	<b>°C</b>	+5 bis +40
Perm. relative humidity	<b>%</b>	80
Protection type acc. to DIN EN 60529		IP 21
Protection class		I (protection earth)
Overvoltage category		II
Contamination level		2
Operation at a terrestrial attitude	<b>m</b>	max. 2000
Dimensions total W x D x H	<b>mm</b>	127 x 149 x 136
Weight without supported load	<b>kg</b>	4.5
Max. supported load incl. attachment	<b>kg</b>	0.5

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**IKA®-Werke GmbH & Co.KG**