

## arium® Water Purification Systems



turning science **into solutions**

# arium<sup>®</sup> Laboratory-grade Water Purification Systems

Our arium<sup>®</sup> laboratory-grade water purification systems inspire with their application-oriented operating designs. They enable you to perform your workflows faster and more reliably, make your daily lab work easier and ensure long-term economical operation. All instruments offer the ultimate flexibility, since they can be perfectly integrated into your laboratory environment and adapted to your requirements.

A choice of more than 70 arium<sup>®</sup> versions is available to meet all your requirements on water quality and to cover any application. The arium<sup>®</sup> bagtanks and the innovative iJust function specially matched to these versions enable significantly more cost-effective water usage and efficient operation of these water purification systems than do conventional units. All arium<sup>®</sup> systems are already certified when delivered. Plus, we offer equipment qualification and maintenance as supplementary services to considerably extend their uses even further.



with Added Value



## Reliable Processes Thanks to the Highest Water Quality

Low detection limits and sensitive analysis devices require laboratory-grade water with consistently high quality. You can always rely on the arium® laboratory-grade water purification systems to provide you with reliable results for your critical and analytical applications. They meet the highest quality requirements and thus ensure the best reproducible results.

## Innovative arium® bagtank System for Time Savings and User Safety

Pure water is stored in the closed arium® bagtank system, which reliably protects purified water from secondary contamination. This ensures consistently high water quality and, therefore, reproducible results.

Unlike conventional tank systems, arium® bagtank technology saves time because the bag can be replaced in less than 5 minutes. At the same time, it increases your users' safety because the bagtank eliminates time-consuming cleaning procedures with hazardous chemicals. The rollers on the arium® bagtank let you move it fast and conveniently to wherever you need to use it.





## Easy Operation Using Touch-activated Functions

You name it, from adjusting basic settings to dispensing water, all functions of the arium® can be controlled by touch screen. Navigate intuitively through the logical user interface – even while wearing gloves. Virtual real-time updates of measured data, flow charts and warning messages are displayed at all times.

Water dispensing can be carried out manually using the slider, the foot switch or via the Favorites function. The latter, in particular, simplifies daily dispensing of recurrent volumes.

## Less Waiting Means More Cost Savings

Consumables, i.e., the filter cartridges, can be quickly and easily changed out. This minimizes maintenance and downtime.

## Optimized Water Usage with iJust

The clever iJust feature automatically optimizes your purified water quality and water usage. Its intelligent software controls a valve on the concentrate drain based on the data measured for  $\text{CaCO}_3$  and  $\text{CO}_2$ .

The advantages:

- Ensures more economical water usage
- Provides the highest purified water quality at all times
- Extends the life of downstream water purification systems

# More Flexibility

The different modules of arium® systems offer the perfect solution for every task in the laboratory. The display is positioned at eye level. You can set up the dispensing location exactly where you need it. Depending on your given space requirements, the system can be integrated as desired at any location within the laboratory.



## Benchtop Unit

Space-saving, compact design; water is dispensed directly at the display level.



## Remote Dispenser with Stand

The remote dispenser features an ergonomic design. Moreover, its height can be adjusted by up to 70 cm (around 30 inches = more than 2 feet). Both of these features enable you to work effortlessly using just one hand. The extended tube guide lets you expand your work area by 3.7 m, or more than 12 ft. This remote dispenser can be used in combination with all benchtop, wall-mounted and built-in units.



arium® pro

## Wall-mounted Unit

The wall-mounted unit saves valuable space on your lab benchtop. The display and dispensing unit are positioned at the bottom to ensure user-friendly, ergonomic operation.



## Built-in Unit

This version saves space on and above your lab bench. You can choose to mount both the display and water dispensing units on a wall or on a multifunctional stand.



## Remote Dispenser with Wall-mounting Plate

This ergonomically designed and easy-to-use remote dispenser is simply mounted on a wall to save space. It is suitable for all benchtop, wall-mounted and built-in units.



## Multifunctional Stand

This convenient stand combines a design that allows unlimited access to all display and dispensing functions and flexibility provided by its height that is adjustable by up to 70 cm (around 30 inches).

# Exactly the Right System – Always

A selection of more than 70 different varieties offers tailor-made solutions without any compromises for all applications.

---

## Features

### **iJust**

Optimized purified water quality and water usage; extends the life of downstream ultrapure water systems

### **arium® bagtank System**

Reliable protection against secondary contamination; easy change-out of the arium® bag saves time; protects users as it eliminates the need for hazardous cleaning chemicals

### **Display with Touch-activated Functions**

Clearly organized user interface featuring intuitive navigation; continuous display of measured data and warning messages

### **High-quality Features**

Integrated feed water conductivity measurement standard on all units; integrated pressure controller, integrated space for depositing utensils, and many more convenient features

### **Maximum System Flexibility**

A variety of hardware configurations; operating and dispensing unit can be flexibly positioned; easy and space-saving integration

### **The Highest Water Quality**

Meets and exceeds ASTM Type 1; excellent retention rates for RNases, DNases and endotoxins, and reduces TOCs; ideal for demanding biological and analytical applications, such as cell cultivation and chromatography

### **SOP Monitoring**

Graphic and acoustic signals for displaying maintenance prompts, alarm messages when limits are exceeded or data falls below minimum limits as well as maintenance interval prompts

### **"Favorites" function**

Quick access to pre-programmed volume dispensing

### **EDI technology**

Consistently high Type 2 water quality; desalted electrochemically

---

## Type of Water

ASTM Type 1

Type 2

Type 3





arium® comfort Series  
Combined pure and ultrapure water systems



arium® pro Series  
Ultrapure water systems;  
any application-specific configuration can be selected



arium® advance Series  
Pure water systems

|  |                          |  |                      |             |
|--|--------------------------|--|----------------------|-------------|
|  | ✓                        |  |                      | ✓           |
|  | ✓                        |  |                      | ✓           |
|  | ✓                        |  | ✓                    | ✓           |
|  | ✓                        |  | ✓                    | ✓           |
|  | ✓                        |  | ✓                    | ✓           |
|  | ✓                        |  | ✓                    |             |
|  | ✓                        |  | ✓                    | ✓           |
|  | ✓                        |  | ✓                    |             |
|  | comfort II               |  |                      | advance EDI |
|  | comfort I and comfort II |  | All arium® pro units |             |
|  | comfort II               |  |                      | advance EDI |
|  | comfort I                |  |                      | advance RO  |



## arium® comfort Series

The arium® comfort lab water systems combine pretreatment and final treatment technologies into one device. Featuring a compact design, they provide ultrapure and pure water of the highest quality. The pure water produced is kept in a closed arium® bagtank. This guarantees optimal storage of the purified water and protects against secondary contamination. The replaceable bag eliminates time-consuming tank cleaning intervals.

The optional UV lamp (185 | 254 nm) built into the ultrapure water loop also prevents microbiological growth and reduces the TOC content to a minimum. The current TOC value is continuously checked by an optionally integrated TOC monitor and displayed for convenient verification.

### Features

#### Safe

TOC content  $\leq 2$  ppb for reproducible results; continuously updated TOC readings can be conveniently viewed on the display

#### Efficient

Optimized water usage ensured by intelligent iJust feature

#### Easy

Glass display with touch-activated functions and intuitive menu navigation

#### Time-saving

The arium® bagtank system provides reliable protection against secondary contamination and eliminates the need for regular tank cleaning

#### Space-saving

Compact design saves valuable work space

#### Fast

"Favorites" function for direct access to pre-programmed volume dispensing

### Critical Application (Type 1)

HPLC

GC-MS, AAS, ICP-MS

Ion chromatography

TOC analysis

SPE (solid phase extraction)

### Standard Applications (Type 2 | Type 3)

Microbiological media and reagents

Histology

ELISA, RIA

Buffers and pH solutions

Feed water for laboratory devices, such as autoclaves, glassware washers, humidifiers, water baths, etc.

### Water Quality

Conductivity for Type 1:  $0.055 \mu\text{S}/\text{cm}$  ( $\approx 18.2 \text{ M}\Omega \times \text{cm}$ )

Typical conductivity for Type 2:  $0.2\text{--}0.07 \mu\text{S}/\text{cm}$  ( $\approx 5\text{--}15 \text{ M}\Omega \times \text{cm}$ )

Typical conductivity for Type 3:  $< 20 \mu\text{S}/\text{cm}$  ( $\approx 0.05 \text{ M}\Omega \times \text{cm}$ )

TOC content for Type 1:  $\leq 2$  ppb

Microorganisms:  $< 1$  CFU/1,000 ml

Particles:  $< 1/\text{ml}$

Type 1 | Type 2 Water

Type 1 | Type 3 Water

## arium® comfort II

The arium® comfort II delivers up to 2 L/min of ASTM Type 1 ultrapure water, as well as 5 L/h or 10 L/h of Type 2 pure water. The additional EDI technology downstream from the RO modules optimizes the system and offers highly efficient ion retention with a quality ranging from 0.2 to 0.07  $\mu\text{S}/\text{cm}$ .

## arium® comfort I

The arium® comfort I provides up to 2 L/min of ASTM Type 1 ultrapure water, as well as 8 L/h or 16 L/h of Type 3 pure water. Feed water is treated with a pretreatment cartridge and downstream RO modules; continuous permeate backflushing effectively prevents deposits and fouling.

### arium® comfort II

- ✓
- ✓
- ✓
- ✓
- ✓

- ✓
- ✓
- ✓
- ✓
- ✓

- ✓
- ✓
- ✓
- ✓
- ✓

### arium® comfort I

- ✓
- ✓
- ✓
- ✓
- ✓

- ✓
- ✓

- ✓
- ✓
- ✓
- ✓
- ✓



Type 1 Water



## arium<sup>®</sup> pro Series

The arium<sup>®</sup> pro ultrapure water systems provide the ultimate flexibility and an excellent cost-benefit ratio because their device configurations are specially tailored to your applications.

All systems meet and exceed the ASTM Type 1 water quality standards and ensure the best reproducible results. Up to 2 L/min of ultrapure water with a conductivity of 0.055  $\mu\text{S}/\text{cm}$  ( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ ) can be dispensed in consistently high quality.

### Features

#### Reliable

TOC content  $\leq 2$  ppb for reproducible results; current TOC readings can be read directly on the display

#### Complete range

Five systems specially customized to meet your specific applications

#### Flexible

Adapts to any laboratory environment thanks to flexible positioning of the system, control unit and dispensing unit

#### Easy

Glass display with touch-activated functions and intuitive menu navigation

#### Fast

"Favorites" function for direct access to pre-programmed volume dispensing



## arium® pro

The arium® pro is a particularly cost-effective system. It focuses on the most important functions and produces ultrapure water in uncompromising quality.

### Standard Applications

- AAS, ICP-MS
- Ion chromatography
- Preparation of reagents
- Photometry

### Water Quality

- Conductivity: 0.055  $\mu\text{S}/\text{cm}$   
( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ )
- TOC content: < 5 ppb
- Microorganisms: < 1 CFU/1,000 ml
- Particles: < 1/ml

## arium® pro DI

The arium® pro DI delivers ultrapure water for standard applications of any kind. The Elemental Kit comprising a set of cartridges reliably removes organic and inorganic components; the TOC content of ultrapure water produced is < 5 ppb.

### Standard Applications

- AAS, ICP-MS
- Ion chromatography
- Preparation of reagents
- Photometry

### Water Quality

- Conductivity: 0.055  $\mu\text{S}/\text{cm}$   
( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ )
- TOC content: < 5 ppb
- Microorganisms: < 1 CFU/1,000 ml
- Particles: < 1/ml

## arium® pro UV

In arium® pro UV, the Analytical Kit cartridges reliably retain organic and inorganic components. The system's integrated UV lamp (185 | 254 nm) prevents microbiological growth and permits TOC values of  $\leq 2$  ppb.

The current TOC value is continuously checked by the optionally integrated TOC monitor and shown on the display.

### Chemical-Analytical Applications

- HPLC
- GC-MS, AAS, ICP-MS
- Ion chromatography
- TOC analysis
- Photometry

### Water Quality

- Conductivity: 0.055  $\mu\text{S}/\text{cm}$   
( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ )
- TOC content:  $\leq 2$  ppb
- Microorganisms: < 1 CFU/1,000 ml
- Particles: < 1/ml



## arium® pro UF

In arium® pro UF, the Biological Kit cartridges reliably remove organic and inorganic components. The system's integrated ultrafilter module ensures that the purified water does not contain any endotoxins, RNA | DNA or DNases and RNases.

The arium® pro UF delivers ultrapure water of consistently high quality and offers ideal conditions to ensure the reliability of critical biological results.

### Chemical-Analytical and Biological Applications:

- AAS, ICP-MS
- Ion chromatography
- Electrophoresis
- Endotoxin analysis
- Immunocytochemistry
- Culture media for cell cultivation
- Production of monoclonal antibodies
- Photometry

### Water Quality

- Conductivity: 0.055  $\mu\text{S}/\text{cm}$  ( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ )
- TOC content: < 5 ppb
- Endotoxins: < 0.001 EU/ml
- RNases: < 0.004 ng/ml
- DNases: < 0.024 pg/ $\mu\text{l}$
- Microorganisms: < 1 CFU/1,000 ml
- Particles: < 1/ml

## arium® pro VF

This high-end arium® pro VF unit delivers constantly high ultrapure water quality. Its integrated UV lamp (185 | 254 nm) and hollow-fiber ultrafilter module not only prevent microbiological growth and reduce the TOC content to  $\leq 2$  ppb, but also remove endotoxins, microorganisms, RNA | DNA and DNases and RNases.

This system is the ideal solution for all critical applications in your laboratory.

### Chemical-Analytical and Biological Applications

- HPLC
- GC-MS, AAS, ICP-MS
- Ion chromatography
- TOC analysis
- PCR
- Electrophoresis
- Endotoxin analysis
- Immunocytochemistry
- Culture media for cell cultivation
- Production of monoclonal antibodies
- Photometry

### Water Quality

- Conductivity: 0.055  $\mu\text{S}/\text{cm}$  ( $\cong 18.2 \text{ M}\Omega \times \text{cm}$ )
- TOC content:  $\leq 2$  ppb
- Endotoxins: < 0.001 EU/ml
- RNases: < 0.004 ng/ml
- DNases: < 0.024 pg/ $\mu\text{l}$
- Microorganisms: < 1 CFU/1,000 ml
- Particles: < 1/ml



## arium® advance Series

The arium® advance EDI and arium® advance RO systems achieve the highest retention rates of ions with an optimal water yield and reliably remove oxidants, heavy metal ions and particles from the feed water. The built-in iJust function also optimizes purified water quality and water usage. Pure water is stored in the innovative, closed arium® bagtank system, which protects this water from secondary contamination.

### Features

#### Efficient

Optimized water usage ensured by the intelligent iJust feature

#### Easy

Glass display with touch-activated functions and intuitive menu navigation

#### Time-saving

The arium® bagtank system provides reliable protection against secondary contamination and eliminates the need for regular tank cleaning

#### Consistently high quality Type 2 water:

State-of-the-art EDI technology

### Standard Applications

---

Microbiological media and reagents

---

Solutions for chemical analysis and synthesis

---

Histology

---

ELISA, RIA

---

Buffers and pH solutions

---

Feed water for various laboratory devices, such as autoclaves, glassware washers, humidifiers, water baths, etc.

### Water Quality

---

Typical conductivity: 0.2–0.07  $\mu\text{S}/\text{cm}$  ( $\cong$  5–15  $\text{M}\Omega \times \text{cm}$ )

---

Typical conductivity: < 20  $\mu\text{S}/\text{cm}$  ( $\cong$  > 0.05  $\text{M}\Omega \times \text{cm}$ )

---

Microorganisms: < 1 CFU/1,000 ml

---

Particles: < 1/ml



Type 2 Water

Type 3 Water

### arium® advance EDI

The arium® advance EDI provides Type 2 pure water in consistently high quality. At a flow rate of 5 L/h or 10 L/h, the system uses the latest EDI technology to ensure reliable and safe removal of impurities in the feed water and reduces the ion content to a minimum.

### arium® advance RO

The arium® advance RO provides Type 3 reverse osmosis water. With a flow rate of 8 L/h, 16 L/h or 24 L/h and automatic RO membrane backflushing, the arium® advance RO is the perfect choice for your general laboratory applications.

#### arium® advance EDI

#### arium® advance RO

|   |   |
|---|---|
| ✓ |   |
| ✓ |   |
| ✓ |   |
| ✓ |   |
| ✓ | ✓ |
| ✓ | ✓ |
|   |   |
| ✓ |   |
|   | ✓ |
| ✓ | ✓ |
| ✓ | ✓ |





# Service for Ultrapure Water Systems

For many applications in the biopharmaceutical industry, at universities and in clinical settings, the availability of pure and ultrapure water of the highest quality is of great importance. Growing demands and more sensitive analytical equipment are raising the bar higher and higher for pure and ultrapure water. Whether you need laboratory-grade water of Type 1, Type 2 or Type 3, Sartorius has the right system for your specific application and the water quality you require.

We invest in the continuous and thorough training of our application and service specialists so that they understand the full range of your applications and can give you well-founded technical advice in accordance with all legal requirements.

## Services Provided

- Installation and startup as well as equipment qualification (IQ | OQ)
- Device configuration based on feed water analysis to achieve the ultimate in ultrapure water quality
- Regular preventive maintenance, including equipment inspection and replacement of consumables
- Extended warranty services

## Your Benefits

- Installation and startup by factory-trained service staff so that you can unleash the full potential of your instruments
- Regular preventive maintenance ensures the reliability of the results and your lab water instrumentation
- Detailed service and calibration records verify compliance with all requirements for documentation

## Sales and Service Contacts

For further contacts, visit [www.sartorius.com](http://www.sartorius.com)

### Europe

#### Germany

Sartorius Lab Instruments  
GmbH & Co. KG  
Weender Landstrasse 94–108  
37075 Goettingen  
Phone +49.551.308.0  
Fax +49.551.308.3289

#### France Et Suisse Romande

Sartorius France  
2, rue Antoine Laurent de Lavoisier  
ZA de la Gaudrée  
91410 Dourdan  
Phone +33.1.70.62.50.00  
Fax +33.1.64.59.76.39

#### Austria

Sartorius Austria GmbH  
Modectenstrasse 22  
1030 Vienna  
Phone +43.1.7965760.0  
Fax +43.1.7965760.24

#### Belgium

Sartorius Belgium N.V.  
Rue Colonel Bourg 105  
1030 Bruxelles  
Phone +32.2.756.06.90  
Fax +32.2.481.84.11

#### Finland Et Baltics

Sartorius Biohit Liquid Handling Oy  
Laippatie 1  
00880 Helsinki  
Phone +358.9.755.951  
Fax +358.9.755.95.200

#### Hungary

Sartorius Hungária Kft.  
Kagyló u. 5.  
2092 Budakeszi  
Phone +3623.457.227  
Fax +3623.457.147

#### Ireland

Sartorius Ireland Ltd.  
Unit 41, The Business Centre  
Stadium Business Park  
Ballycolin Road  
Dublin 11  
Phone +353.1.8089050  
Fax +353.1.8089388

#### Italy

Sartorius Italy S.r.l.  
Viale A. Casati, 4  
20835 Muggiò (MB)  
Phone +39.039.4659.1  
Fax +39.039.4659.88

#### Netherlands

Sartorius Netherlands B.V.  
Phone +31.30.60.53.001  
Fax +31.30.60.52.917  
[info.netherlands@sartorius.com](mailto:info.netherlands@sartorius.com)

#### Poland

Sartorius Poland sp.z o.o.  
ul. Wrzesinska 70  
62-025 Kostrzyn  
Phone +48.61.6473830  
Fax +48.61.6473839

#### Russian Federation

LLC "Sartorius ICR" and LLC "Biohit"  
Uralskaya str. 4, Lit. B  
199155, Saint-Petersburg  
Phone +7.812.327.5.327  
Fax +7.812.327.5.323

#### Spain Et Portugal

Sartorius Spain, S.A.  
Avda. de la Industria, 32  
Edificio PAYMA  
28108 Alcobendas (Madrid)  
Phone Spain +34.902.123.367  
Phone Portugal +351.800.855.800  
Fax Spain +34.91.358.96.23  
Fax Portugal +351.800.855.799

#### Switzerland

Sartorius Mechatronics Switzerland AG  
Ringstrasse 24a  
8317 Tagelswangen (ZH)  
Phone +41.44.746.50.00  
Fax +41.44.746.50.50

#### U.K.

Sartorius UK Ltd.  
Longmead Business Centre  
Blenheim Road, Epsom  
Surrey KT19 9QQ  
Phone +44.1372.737159  
Fax +44.1372.726171

#### Ukraine

LLC "Biohit"  
Post Box 440 "B"  
01001 Kiev, Ukraine  
Phone +380.44.411.4918  
Fax +380.50.623.3162

### Americas

#### USA

Sartorius Corporation  
5 Orville Drive, Suite 200  
Bohemia, NY 11716  
Phone +1.631.254.4249  
Toll-free +1.800.635.2906  
Fax +1.631.254.4253

#### Argentina

Sartorius Argentina S.A.  
Int. A. Avalos 4251  
B1605ECS Munro  
Buenos Aires  
Phone +54.11.4721.0505  
Fax +54.11.4762.2333

#### Brazil

Sartorius do Brasil Ltda  
Avenida Senador Vergueiro 2962  
São Bernardo do Campo  
CEP 09600-000 - SP- Brasil  
Phone +55.11.4362.8900  
Fax +55.11.4362.8901

#### Canada

Sartorius Canada Inc.  
2179 Dunwin Drive #4  
Mississauga, ON L5L 1X2  
Phone +1.905.569.7977  
Toll-Free +1.800.668.4234  
Fax +1.905.569.7021

#### Mexico

Sartorius de México S.A. de C.V.  
Circuito Circunvalación Poniente  
No. 149  
Ciudad Satélite  
53100, Estado de México  
México  
Phone +52.5555.62.1102  
Fax +52.5555.62.2942

### Asia | Pacific

#### Australia

Sartorius Australia Pty. Ltd.  
Unit 5, 7-11 Rodeo Drive  
Dandenong South Vic 3175  
Phone +61.3.8762.1800  
Fax +61.3.8762.1828

#### China

Sartorius Scientific  
Instruments (Beijing) Co., Ltd.  
33 Yu An Road, Airport Industrial Park  
Zone B,  
Shunyi District, Beijing 101300,  
P.R.China  
Phone +86.10.8042.6300  
Fax +86.10.8042.6486

#### Hong Kong

Sartorius Hong Kong Ltd.  
Unit 1012, Lu Plaza  
2 Wing Yip Street  
Kwun Tong  
Kowloon, Hong Kong  
Phone +852.2774.2678  
Fax +852.2766.3526

#### India

Sartorius Weighing India Pvt. Ltd.  
#69/2-69/3, NH 48, Jakkasandra,  
Nelamangala Tq  
562 123 Bangalore, India  
Phone +91.80.4350.5250  
Fax +91.80.4350.5253

#### Japan

Sartorius Japan K.K.  
4th Fl., Daiwa Shinagawa North Bldg.  
8-11, Kita-Shinagawa 1-chome  
Shinagawa-ku, Tokyo, 140-0001 Japan  
Phone +81.3.3740.5408  
Fax +81.3.3740.5406

#### Malaysia

Sartorius Malaysia Sdn. Bhd  
Lot L3-E-3B, Enterprise 4  
Technology Park Malaysia  
Bukit Jalil  
57000 Kuala Lumpur, Malaysia  
Phone +60.3.8996.0622  
Fax +60.3.8996.0755

#### Singapore

Sartorius Singapore Pte. Ltd  
1 Science Park Road,  
The Capricorn, #05-08A,  
Singapore Science Park II  
Singapore 117528  
Phone +65.6872.3966  
Fax +65.6778.2494

#### South Korea

Sartorius Korea Ltd.  
8th Floor, Solid Space B/D,  
PanGyoYeok-Ro 220, Bundang-Gu  
SeongNam-Si, GyeongGi-Do, 463-400  
Phone +82.31.622.5700  
Fax +82.31.622.5799

#### Thailand

Sartorius (Thailand) Co. Ltd.  
129 Rama 9 Road,  
Huaykwang  
Bangkok 10310  
Phone +66.2643.8361-6  
Fax +66.2643.8367



◀ [www.sartorius.com](http://www.sartorius.com)