

Product Name CHEMGENE MEDLAB Multi-Surface Disinfectant Wipes – 96 Pack

Product Overview CHEMGENE MEDLAB multi-surface disinfectant wipes are designed for hard surface applications in laboratory, critical services, and forensics settings.

Intended to provide convenience of application without compromising antimicrobial efficacy, CHEMGENE wipes work to decontaminate surfaces and equipment in just 60 seconds, removing dirt and debris and denaturing DNA/RNA.

CHEMGENE MEDLAB multi-surface disinfectant wipes are manufactured to strict ISO 9001 quality standards in the UK, and are compliant with REACH, CLP and BPR regulations in the UK. The plastic-free wipe substrate is biodegradable within 3 months.


For more information, technical documentation and training materials, visit byotrol.com. Use biocides safely. Always read the label and product information before use.

Microbiological Testing CHEMGENE MEDLAB multi-surface wipes have been tested to be effective under the following microbiological test conditions: -

EN 13727 (bactericidal)
EN 16615 (bactericidal – disinfectant wipe test)
EN 13623 (Legionella)
EN 13624 (yeasticidal)
EN 16615 (yeasticidal – disinfectant wipe test)
EN 13624 (fungicidal)
EN 14476 (virucidal against enveloped viruses)

Added Benefits

- Denatures DNA/RNA in 5 minutes
- Biodegradable formulation
- Biodegradable, plastic-free wipe substrate
- Manufactured in the UK
- Non-corrosive
- Alcohol-free
- No rinsing required
- Compatible with a wide range of materials
- Compliant with REACH, CLP and biocidal regulations in the UK and EU
- 2 year shelf life

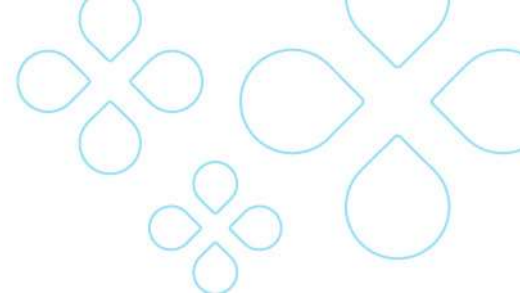

Directions for Use

Suitable for use on hard, non-porous surfaces. Remove any visible dirt or soiling before disinfecting. Lift seal and remove one wipe from pack. Place palm of hand on top of the wipe. Apply the wipe to the surface in an 'S' pattern, ensuring full coverage. Each wipe is suitable for disinfecting 1m². Avoid wiping the same area twice and use a new wipe for each new surface, to avoid cross-contamination. Leave surfaces to dry naturally, no need to rinse. Reseal pack after use to avoid wipes drying out. Test an inconspicuous area before use if unsure. Not to be used to disinfect medical devices.

Dispose of used wipes in general waste or clinical waste as appropriate. Wipes are biodegradable within 3 months.

Hazard Class

Non-hazardous

PPE Requirements

None

Byotrol Technology Ltd

Address:
Building 303 (Ashton), Thornton Science Park,
Pool Lane, Chester, CH2 4NU, UK

Email:
info@byotrol.com

Web:
www.byotrol.com

Telephone:
+44 (0) 1925 742 000

LinkedIn:
 [Byotrol Plc](#)