

The following

POLYSAFE®

Terry knitted para-aramid nitrile lined glove series are manufactured by
Bennett Safetywear Ltd, (BSL)

STOCKCODE DESCRIPTION SIZE

1515900011	FTK/NT/25	9
1515950031	FTK/NT/35	9
1515970021	FTK/NT/45	9



Gloves are made to BS EN ISO 21420:2020 sizes.

CE 0598 UK CA 0120

STANDARD EN 388:2016 + A1:2018

	Level
Abrasion Resistance	4
Cut Resistance	5
Tear Resistance	4
Puncture Resistance	4
Cut Resistance (TDM)	E



EN 407:2020

Limited Flame Spread	4
Contact Heat	3
Convective Heat	4
Radiant Heat	3
Small Drops of molten Metal	X
Large Quantities of Molten Metal	X



GLOVE DESCRIPTION

100% Para-aramid outer terry knitted glove lined with nitrile dipped cotton inner. Gloves offer excellent cut resistance and excellent thermal insulation when handling hot objects.

MARKING

Each packet shall be labelled / contain the appropriate marking with regards to the amount, type, size, manufacturer, performance levels and UKCA &

BENNETT SAFETYWEAR LTD

7-11 MERSEY ROAD, CROSBY, LIVERPOOL, UK

TD NO : 1024

CE mark. Labels on the gloves will contain similar information. This information will be provided in the minimum resale quantity.

COMPLEX SAFETY CATERGORY

Gloves are designed to protect hands in the working environment in accordance with EN388:2016 + A1:2018, EN ISO 21420:2020 & EN407:2020. When selecting a glove based on risk analysis it should be understood that the protection is limited to the risk level and standards mentioned above.

The type examination was carried out by ITS Testing Services, Centre Court, Meridian Business Park, Leicester, LE19 1WD, UK. Their test house number is 0362

None of the materials or processes used in the manufacture of these products is known to be harmful to the wearer.

WARNINGS:

Test results apply to the gloves in the as received condition but do not extend to any leather cuff. Result may differ if cleaned. Do not use near moving machinery due to entanglement hazard. Overall classification may not reflect the performance of only the outermost layer. These gloves may not be suitable for protection against sharply pointed objects such as hypodermic needles.

CARE / MAINTENANCE

Cleaning and disinfection is not intended for these gloves.

Both new and used gloves should be thoroughly inspected before and after use, especially after cleaning treatment, and before being worn to ensure no damage is present. Damaged gloves should not be worn and should be disposed of. Wet or contaminated gloves should not be used. Dirty gloves may not provide the same level of protection as that shown. Gloves should not be left in a contaminated condition if re-use is intended, in which case gloves should be cleaned as far as possible, provided no serious hazards exist, before removal from the hands. Excess contaminant should first be removed, e.g. loose dirt can be brushed off with a soft brush. The gloves may be de-contaminated with mild detergent solution, then rinsed with clean water and allowed to dry naturally, ideally with some air movement. When the contaminant is not removable or presents a potential hazard it is advisable to ease left and right hand gloves off alternatively using the gloved hand so that the gloves are removed without the contaminant contacting the bare hands. Washing is not recommended.

Do not wring. Do not tumble dry. Do not use bleach. Gloves may be rinsed in water and allowed to drip dry in ambient temperatures. Reshape whilst still damp.

Warning: 

PACKING AND STORAGE

Can be used under normal climatic conditions.
Gloves shall be wrapped in polymeric packaging.

Gloves can be used up to five years after date of supply if stored correctly. Store in original packaging. Gloves should be ideally stored in a cool dry place away from direct sunlight at ambient temperatures (between 5°C and 25°C) in a dry well ventilated area in original packaging to maintain the optimum properties of the glove. The gloves are packed in bundles, along with this leaflet. This bundle is suitable for transportation and storage.

OBSOLESCENCE

Stored correctly, the gloves' physical properties will not change for up to five years.

GENERAL

The quality systems used to manufacture and supply the gloves are in compliance with ISO 9001:2015.

PLEASE NOTE

The information contained here is intended to assist the wearer in the selection of personal protective equipment. The result of the laboratory tests should help with correct glove selection; however, it should be understood that the actual conditions of use cannot be directly simulated. It is therefore the responsibility of the end user and not the manufacturer to determine the glove suitable for the intended use.

DECLARATION

Bennett Safetywear Ltd declares that the new PPE as described in the TD 1024 technical specification for POLYSAFE FTK/NT GLOVES are in conformity with Regulation (EU) 2016/425 and as amended to apply in GB and with the national standard transposing harmonised standard No : BS EN 407:2020, BS EN 388:2016 + A1:2018, BS EN ISO 21420:2020, and is identical to the PPE which is subject of UKCA certificate LECF00384920 issued by: I.T.S. TESTING SERVICES, CENTRE COURT, MERIDAN BUSINESS PARK, LEICESTER, U.K., LE19 1 WD, Approved Body Number: AB0362 who performed the type-examination and is subject to the procedure set out in the legislation Regulation (EU) 2016/425 and as amended to apply in GB Module D under the supervision of the approved body SGS United Kingdom, Rossmore Business Park, Ellesmere Port, South Wirral, Cheshire, CH65 3EN, United Kingdom. Their approved body number is 0120 and is also the subject of the EU certificate. No: and identical to the PPE which is the subject of the EU certificate. No: LECF00372385, issued by: I.T.S. TESTING SERVICES, CENTRE COURT, MERIDAN BUSINESS PARK, LEICESTER, U.K., LE19 1 WD, Notified Body Number: 0362 by INTERTEK Italia S.p.A. Via Guido Miglioli 2/A 20063 Cernusco sul Naviglio - Milano (MI) Italy NB 257 now owns this certificate., and is subject to the procedures under the supervision of the notified body SGS Fimko Oy, P.O. Box 30 (Särkiniementie 3), 00211 HELSINKI, Finland 0598. Their notified body number is 0598. Done at : Bennett Safetywear Ltd, 7-11 MERSEY ROAD, CROSBY, LIVERPOOL, UK

Further copies of this information and declarations of conformity are available on request from:

Bennett Safetywear Ltd

+44 (0)151 924 3996

Web : <http://bennettsafetywear.co.uk>